

ULVAC

Instruction Manual
for
Direct-Drive Oil Sealed Rotary Vacuum Pump

Model

G-101D

G-101S

Before using the product, be sure to read this manual.
Keep this manual in a place where it can be referred to at any time
and look after it carefully.
The contents of this instruction manual are subject to change
without prior notice due to improvements in performance and the
functions of the product.

ULVAC KIKO,Inc

0. Introduction

0.1 Before using the vacuum pump

Thank you for purchasing our vacuum pump (hereinafter called “pump”). When you have received the pump, check that the delivered pump is as per your order and that it has not been damaged in transportation, etc.



Warning

In order to use the pump for as long as possible, read this instruction manual thoroughly before performing installation, operation, inspection and maintenance, and sufficiently understand the cautions for safety, the specifications and operation methods of the pump.



Note

No part of this instruction manual may be copied for use by a third party without our permission.

0.2 Safety symbols

In this instruction manual and on warning labels attached to the pump, the following symbols are used so that matters which must be strictly adhered to can be readily understood.

These symbols are divided as shown below.

Danger

When mishandled, there is an imminent danger of the operator suffering a fatal accident or serious injury.

Warning

When mishandled, there is a possibility of the operator suffering a fatal accident or serious injury.

Caution

When mishandled, there is a possibility of the operator suffering an injury (light or medium injury) or of damage occurring to property.

Note

When mishandled, there is a possibility of the pump being damaged or malfunctioning.

0.3 Cautions for safety

Danger

When toxic or flammable gases are exhausted from the pump, they may leak not only from the pump outlet but also from the pump unit itself. Take proper measures suitable for the type of gas.

Danger

After the pump has been used for exhausting toxic gasses, not only the pump itself but also vacuum pump oil (hereinafter called “pump oil”) get toxic. Keep this in mind when performing the maintenance.

Warning

Never allow people other than repair engineers to disassemble or repair the pump. Failure to do so may result in ignition or malfunction, leading to injury or electric shock.

Warning

Before performing inspection or repair, always turn off the power switch. Failure to do so may result in electric shock or the unexpected start of the pump, leading to injury.

Warning

Connect the earth wire correctly. It is recommended that a dedicated earth leakage breaker should be installed. If the earth wire is not connected, there is a possibility of electric shock occurring in the case of a malfunction or electrical leakage.

Warning

There is a risk of explosion. Never block the outlet or operate the pump with equipment mounted at the outlet side which blocks the passage of gas. Otherwise, the internal pump pressure increases causing the pump to explode, the oil level gauge to protrude or the motor to be overloaded. This pump is not resistant to pressure. The internal pump pressure is limited to 0.03 MPa (gauge pressure).

 **Warning**

Do not use the pump in an explosive atmosphere. Failure to do so will result in injury or fire.

 **Caution**

Do not insert fingers or objects into the opening of the motor. Doing so may result in electric shock, injury, or fire.

 **Caution**

Never touch the rotating section of the motor, shaft or coupling while the pump is in operation. Failure to do so will result in injury.

 **Caution**

Never place combustible materials around the motor or pump. There is a risk of fire.
Also, do not place objects which block ventilation around the motor. Abnormal heat generation may result in burns or fire.

 **Caution**

Do not touch the motor while the pump is in operation or when the pump is still hot immediately after it stops. Touching it will result in burns.

 **Caution**

Arrange wires correctly in accordance with the “Electrical Equipment Technical Standard” and “Wiring Regulations.” Incorrect wiring may result in fire.

 **Caution**

If the pump ceases operation or malfunctions, turn off the power switch immediately to prevent accidents, and ask the company from which you purchased the pump or the manufacturer for inspection and repair.

 **Note**

Do not operate the pump without adding pump oil. If it is operated in an oil-less condition, the pump will be damaged.

0.4 Acceptance and storage of the pump

0.4.1 Acceptance of the pump

Although the pump is delivered with great care, check the following after unpacking.

- ① The delivered pump is in accordance with your request.
- ② The specified accessories (enough pump oil to use the pump once; optional equipment) have been provided.
- ③ No parts have been damaged in transportation.
- ④ Neither screws nor nuts have become loose nor were lost in transportation.

If there are any problems, contact the company from which you purchased the pump or the sales department of the manufacturer.

0.4.2 Environmental conditions for storage, installation and operation

Since this pump is precisely engineered, ensure that the following conditions be satisfied during storage, installation and operation.

- ① Ambient temperature, relative humidity: 7°C ~ 40°C, 85% RH or less
- ② Height above sea level during storage and installation: 1,000 m or less
- ③ Other conditions for storage and operation
 - a) Free from corrosive and explosive gases
 - b) No condensation
 - c) Dust-free environment
 - d) Indoors
 - e) Do not place pumps on top of each other or place a pump on its side.
 - f) Not subject to direct sunlight
 - g) Far from heat sources
 - h) When you keep it for a long period of time, put pump oil into a pump and seal a inlet port with a cap.
 - i) Don't keep it, where moisture is attracted.

Caution

Since the pump weighs as much as 20 kg, do not lift or transport it by yourself. Doing so may cause an injury. Please wear safety shoes at the time of work. Perform such work by two people as shown in 3.1 "Installation." and wear safety shoes.

Note

Do not subject the pump to shocks or place the pump on its side. Doing so may damage the pump.

Note

For indoors use only.

 **Note**

When you keep it, without using a pump for a long period of time, please put pump oil into a pump and seal a inlet port. Oil is not put into a pump, but if it is kept where a inlet port is opened wide, water absorption may expand vanes of a pump and a pump may stop rotating.

 **Note**

Please do not leave it, where moisture is sucked. If it is left with moisture sucked, since water absorption expansion of vanes and the corrosion of pump parts will be promoted, a pump may stop rotating.

0.5 Protective device

The pump is provided with a single-phase motor.

An overload protector (manually reset thermal protector) is incorporated.

The use of another protective device (such as an earth leakage breaker) in addition to the overload protector is recommended.

 **Note**

Use the pump only at the rated voltage. Use at other than the rated voltage will interfere with correct operation of the overload protector, and result in the motor burning out, or fire.

Contents

0. Introduction	01
0.1 Before using the vacuum pump	01
0.2 Safety symbols	02
0.3 Cautions for safety	03
0.4 Acceptance and storage of the pump	05
0.4.1 Acceptance of the pump	05
0.4.2 Environmental conditions for storage, installation and operation	05
0.5 Protective device	06
1. For Safe Operation	1
1.1 Hazards peculiar to the pump and safety measures	1
1.1.1  Danger Leakage of hazardous gases and substances	1
1.1.2  Warning Electric shock	1
1.1.3  Warning Explosion	2
1.1.4  Caution High temperature	2
1.2 Safety Data Sheet (SDS)	2
2. Outline of the Pump	3
2.1 Specification	3
2.2 Dimensional drawing	5
3. Installation	6
3.1 Installation	6
3.2 Lubrication	6
3.3 Vacuum piping	8
3.4 Electric wiring	9
3.5 Fluctuation in the power voltage and frequency	11
4. Operation	13
4.1 Cautions for operation	13
4.2 Start of operation	14
4.3 Stopping the operation	14
4.4 Backflow preventer (G-101S)	15
4.5 Thermal protector	15
4.6 Gas ballast valve (G-101D)	15
4.7 Operation in cold climates	16
4.7.1 Operational method	16
4.7.2 Important points	16
4.8 An inlet port, exhaust port replacement (at option)	17
4.8.1 The switching method of the inlet port	17
4.8.2 The switching method of the exhaust port	17
4.9 Installation of the oil mist trap (Option)	17
4.10 Restriction on operation when the oil mist trap is installed	17

5. Pump Performance	18
5.1 Ultimate pressure	18
5.2 Pumping speed	18
5.3 Power requirement	18
6. Maintenance, Inspection and Repair	20
6.1 Maintenance	20
6.2 Periodic inspection	20
6.3 Replacement of the pump oil	22
6.4 Replacement of the coupling	24
6.5 Trouble check list	25
7. Disposal	27
8. Maintenance parts	28
8.1 Maintenance parts list	28
8.2 Disassembly drawing	30

Warranty

Safety Data Sheet (SDS)

Pump Usage Check Sheet (Use this sheet for requesting an overhaul.)

Contact address of sales and service departments

Figures and Tables

Fig. 1	Dimensional drawing of G-101D oil sealed rotary vacuum pump	5
Fig. 2	Dimensional drawing of G-101S oil sealed rotary vacuum pump	5
Fig. 3	Transportation method of the oil sealed rotary vacuum pump	6
Fig. 4	Lubrication of the oil sealed rotary vacuum pump	7
Fig. 5	Basic piping diagram to the vacuum chamber	8
Fig. 6	Electric wiring diagram	9
Fig. 7	Change region of the voltage and frequency	11
Fig. 8	Pumping speed curve	19
Fig. 9	Replacement of the coupling	24
Fig. 10	Disassembly drawing of G-101D oil sealed rotary vacuum pump	30
Fig. 11	Disassembly drawing of G-101S oil sealed rotary vacuum pump	31
Table 1	Pump specification	3
Table 2	Motor specification	4
Table 3	Characteristics of the thermal protector	15
Table 4	Periodic inspection table	21
Table 5	Trouble check list	25
Table 6	G-101D Maintenance parts list	28
Table 7	G-101S Maintenance parts list	29
Attached table:	Safety Data Sheet (SDS)	

1. For Safe Operation

1.1 Hazards peculiar to the pump and safety measures

Before operating or inspecting the pump, read this section carefully to fully understand potential hazards and prevention methods.

1.1.1 Danger Leakage of hazardous gases and substances

Cause	Prevention method and measures
Leakage of toxic and flammable gases	⇒ Dilute hazardous gases to a safe concentration before they enter the pump inlet.
Injury due to touching toxic pump oil in the pump or harmful substances attached to the pump during inspection or disposal	⇒ ① Wear protective equipment suitable for toxic substances when carrying out inspection. ② Before overhauling and disposing of the pump, ask a waste disposal specialist to make it safe. ③ Ask an authorized waste disposal specialist to carry out disposal.

1.1.2 Warning Electric shock

Cause	Prevention method and measures
The energized part of the motor was touched.	⇒ ① When connecting electric wires, always turn off the power and be sure to connect the earth wire. ② When inspecting and transferring the pump, always turn off the power. ③ Never insert hands, fingers, or thin objects through the motor opening.

1.1.3 Warning Explosion

Cause	Prevention method and measures
The pressure in the pump increased causing the pump to explode.	⇒ The maximum internal pump pressure is 0.03 MPa (gauge pressure). Measure the pressure at the outlet side and, if the pressure is 0.03 MPa or more (gauge pressure), remove objects which block the passage of gas from the outlet side. When an oil mist trap is adopted, replace or clean it so that it will not block the passage of gas.

1.1.4 Caution High temperature

Cause	Prevention method and measures
High temperatures caused burns.	⇒ ① The pump reaches a high temperature during operation. pump body → 70 ~ 80 °C motor → 70 ~ 80 °C ② Since the surface temperature is hot, touching the surface accidentally may result in burns. Never touch the pump during operation. When carrying out inspection, wait until the pump has cooled down completely after it stops.

1.2 Safety Data Sheet (SDS)

The attached “Safety Data Sheet (SDS)” shows chemical materials which may be used or touched when operating the pump. Read the SDS carefully in order to understand the harmful properties of these materials.

Contact us before using chemical materials (vacuum pump oil) other than those mentioned in this instruction manual.

Caution

SDS is submitted as reference information to ensure safe handling of hazardous and harmful materials. Personnel handling the pump oil should be aware that proper measures must be taken depending on the conditions of use as their responsibility. Keep in mind that the SDS itself is not a warranty for safety.

2. Outline of the Pump

2.1 Specification

This oil sealed rotary vacuum pump is a rotary vane pump (hereinafter called Gaede type pump) in which the pump is directly driven by the motor. Since the pump is small, light, and quite simply constructed, it is easily maintained and repaired.

Table 1 Pump specification

Model		G-101D	G-101S
Type		Sliding vane 2stage	Sliding vane 1stage
Pumping speed (L/min)	50Hz	100	
	60Hz	120	
Ultimate pressure (Pa)	G.V. Close	6.7×10^{-1}	9.3
	G.V. Open	6.7	-
Oil	Standard oil	SMR-100	
	Oil amount (mL)	800	1,200
Inlet port diameter		$\phi 27$ (Inlet port of 40□Flange can be installed in pump)	
Exhaust port diameter		Mushroom type (Exhaust port of G1size screw can be installed in pump)	
Weight (Kg)	Single phase	23	22

Note 1: The ultimate pressure values in the above table are indicated by a Pirani gauge.

Note 2: Vacuum pump oils have different steam pressures, viscosities, and oil properties depending on the type. Always use the oil sealed rotary vacuum pump oil specified by us. The use of other oils will affect the pump's performance.

Specified oil: SMR-100

Note 3: "G.V." is an abbreviation for gas ballast valve.

Table 2 Motor specification

Model			G-101D	G-101S
Motor	W(Poles)		400(4)	
	Single phase	V	100/100(50/60Hz)	
			200/200(50/60Hz)	
Full load current (A)	Single phase	100/100V	6.4/5.9(50/60Hz)	
		200/200V	3.2/2.8(50/60Hz)	
Motor speed (r/min)	Single phase	100/100V	1,425/1,710(50/60Hz)	
		200/200V	1,450/1,745(50/60Hz)	



Caution


Do not touch the motor while the pump is in operation or when the pump is still hot immediately after it stops. Touching it will result in burns.



Note

Use the pump only at the rated voltage. Use at other than the rated voltage will interfere with correct operation of the overload protector, and result in the motor burning out, or fire.


2.2 Dimensional drawing

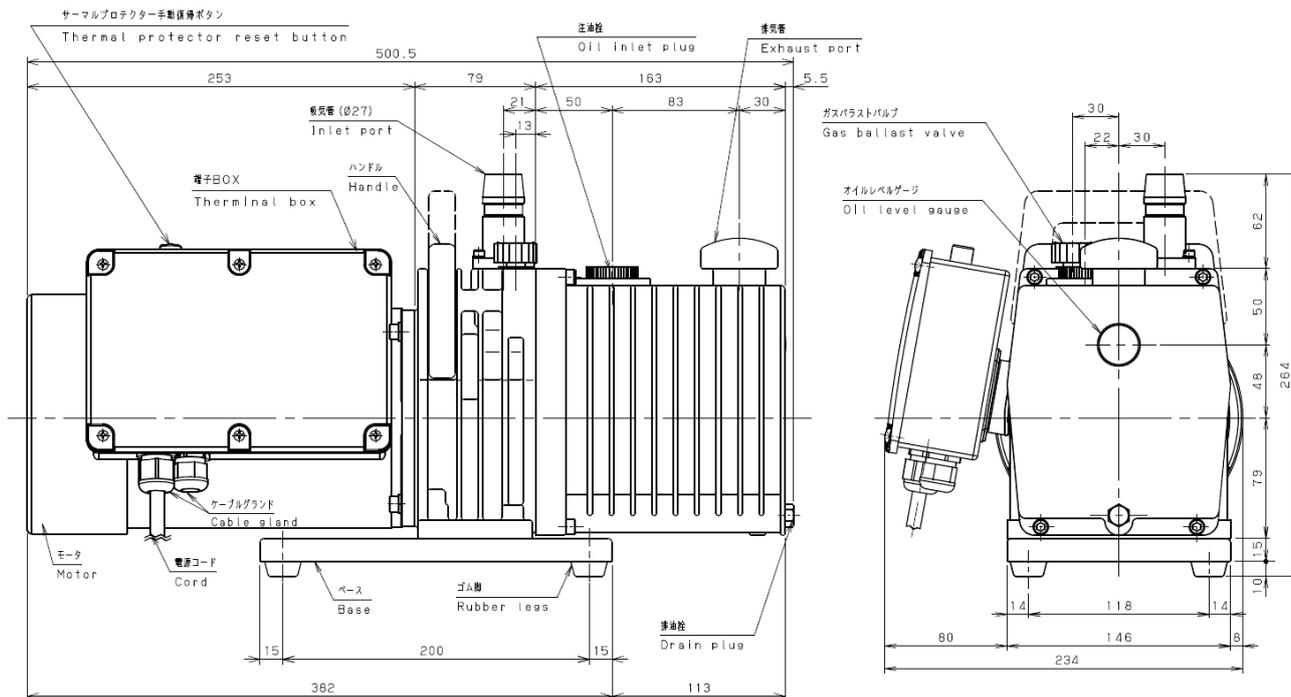

Motor ; Single phase, 50/60Hz, 400W, 4Poles, Capacitor start & run,
fully-closed external fan

Weight ; 23kg

Fig. 1 Dimensional drawing of G-101D oil sealed rotary vacuum pump

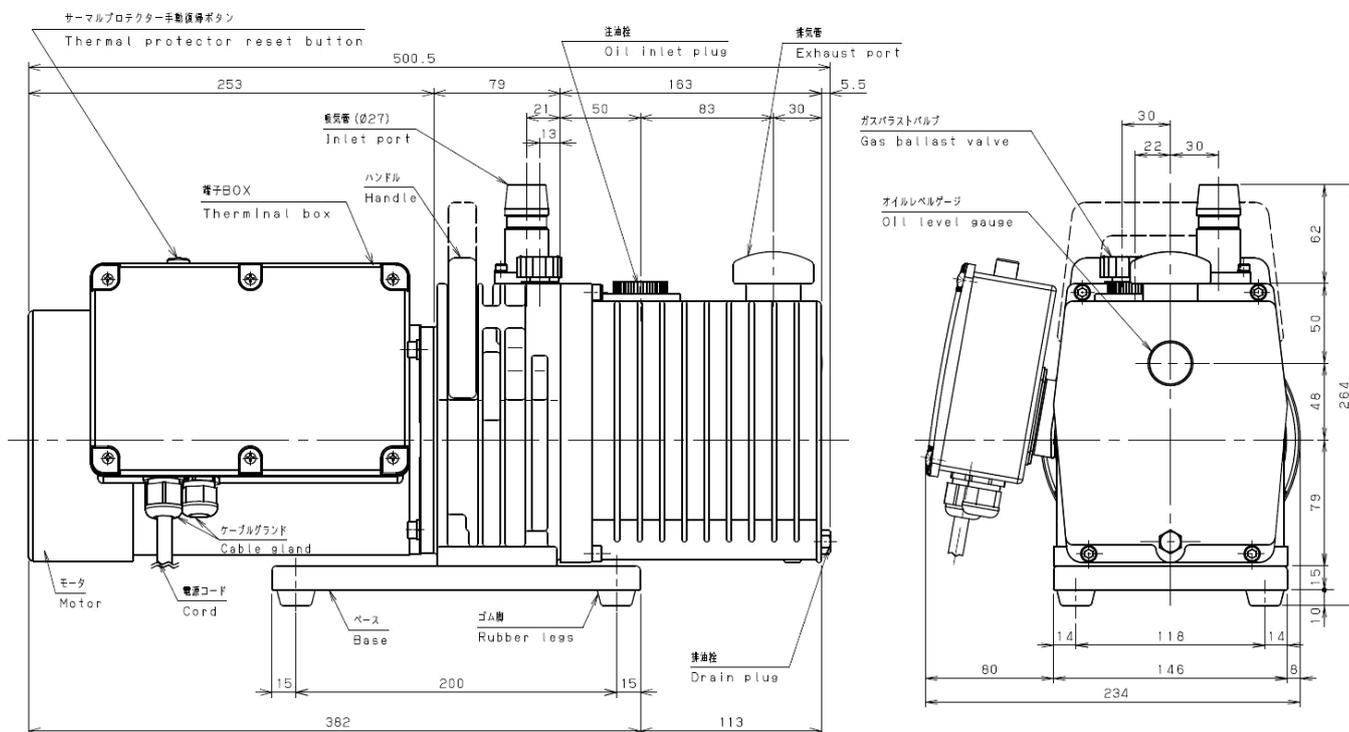

Motor ; Single phase, 50/60Hz, 400W, 4Poles, Capacitor start & run,
fully-closed external fan

Weight ; 22kg

Fig. 2 Dimensional drawing of G-101S oil sealed rotary vacuum pump

3. Installation

3.1 Installation

The pump should be installed on a level surface in a location with minimal dust, dirt and humidity and be arranged with consideration given to ease of installation, removal, inspection and cleaning.

Particular attention should be paid to the ambient temperature when building the pump into equipment. Use a rubber vibration isolator to separate the pump from other equipment and to isolate the pump from the vibrations of other equipment. See “0.4.2 Environmental conditions for installation, storage and operation” for details.



Fig. 3 Transportation method of the oil sealed rotary vacuum pump

Caution

Since the pump weighs as much as 20 kg, do not lift or transport it by yourself. Doing so may cause an injury. Always carry out work with two people as shown in Fig. 3.

Note

If the pump is operated whilst it is tilted, placed on its side or upside-down, the pump will be damaged. Install the pump level with the inlet facing up as shown in Fig. 1, Fig. 2.

3.2 Lubrication

Remove the lubrication plug from the lubrication port, and add the pump oil which has been delivered together with the pump or the pump oil specified by us (SMR-100) up to the range marked with the red line on the oil level gauge. When making the first lubrication, add oil near to the upper oil level limit shown on the oil level gauge. After lubrication, mount the lubrication plug to the pump (see Fig. 4).

Always keep the oil level of the pump within the oil limit range shown on the oil level gauge during operation. If the amount of oil is incorrect, the performance of the pump will deteriorate resulting in the malfunctioning of the pump. When the amount of oil has reduced and the oil level has reached an area below the lower red line which shows the lower limit on the oil level gauge such that the level cannot be seen, the ultimate pressure increases and exhausting sound may not cease.

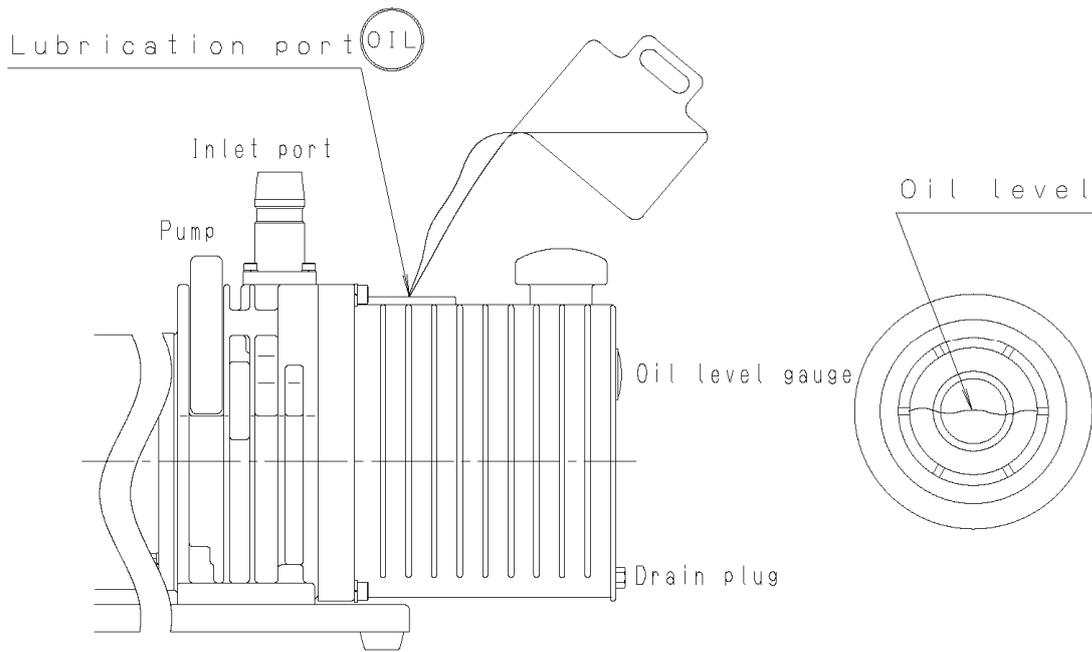


Fig. 4 Lubrication of the oil sealed rotary vacuum pump

⚠ Caution

- ① Wear protective equipment such as rubber gloves and safety goggles.
- ② Be sure to read the attached "Safety Data Sheet" before adding oil. If the oil accidentally comes into contact with your hands or enters your eyes, take proper measures in accordance with the section "First-aid treatment" shown in "Safety Data Sheet."

⚠ Note

Use only oils specified by us. If other oils are used, the pump performance will deteriorate or its life will be shortened.

3.3 Vacuum piping

- (1) Before connecting the pipe to the pump, clean the inner walls of the vacuum chamber, piping and vacuum valve to completely eliminate moisture, fine particles, dust, dirt and rust.

⚠ Note

If fine particles, dust or dirt, etc are evacuated, the pump may malfunction. If moisture is evacuated, not only does the ultimate pressure increase but also the inside of the pump becomes rusty causing the pump to malfunction.

- (2) Mount vacuum valve (A) and leak valve (B) between the vacuum chamber and pump as shown in Fig. 5

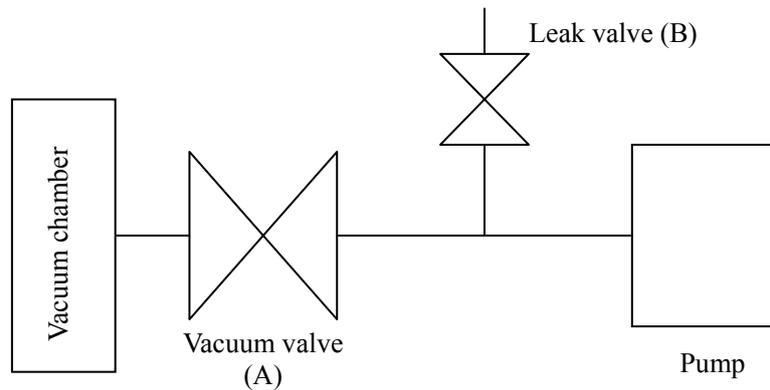


Fig. 5 Basic piping diagram to the vacuum chamber

- (3) Vacuum tubing between inlet port and another port is provided with vacuum rubber hose.

⚠ Note

The wire mesh in the inlet port has been adopted to prevent foreign matter from entering the pump. Do not remove the wire mesh.

3.4 Electric wiring

- (1) This pump has performed beforehand electric wiring by the side of a pump.
When wiring, open the terminal box of the motor and connect the wires as shown in Fig. 6.
- (2) Confirm that use power supply voltage accords with a working voltage of the motor.
- (3) The pump rotates in the clockwise direction as seen from the front of the pump (level gauge side).
- (4) Fault load protection equipment (manual return type thermal protector) is built in this motor.
- (5) The use of another protective device (such as an earth leakage breaker) in addition to the overload protector is recommended.
- (6) This motor can pick up the operation of the thermal protector as signaling.

A lead wire for thermal sign takeoff is connected to a terminal block in the terminal box.

When I take out thermal sign, I open terminal box of the motor, and please connect it like fig. 6.

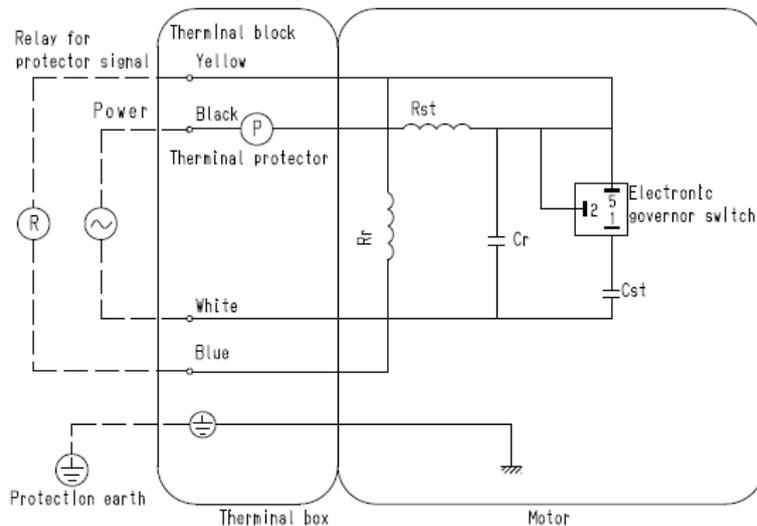


Fig. 6 Electric wiring diagram

Warning

A start mode of this motor is condenser start / operating method. The switching of the start / operation reaches by electronic governor switch. When I apply voltage of the inverter control because I use an electron element for electronic governor switch, malfunction of the electronic governor switch occurs and causes the motor stumbling-block. Please never apply voltage of the inverter control.

 **Warning**

Before connecting wires, turn off the power switch. Never perform wiring with the power supplied as an electric shock will occur. Connect the earth wire correctly. Failure to do so may result in electric shock if a failure or earth leakage occurs. Installation of a dedicated earth leakage breaker is also recommended.

 **Note**

Perform electric wiring correctly in accordance with the “Electric Equipment Technical Standard” and “Internal Wiring Regulation.” Incorrect wiring will result in fire.

 **Warning**

If the pump is directly (and permanently) connected to the host equipment then the end-user has to provide adequate disconnection device.

 **Caution**

GROUNDING INSTRUCTIONS

In the event of an electrical short circuit, grounding reduces the risk of electric shock by providing an escape wire for the electric current. This pump is equipped with a cord having a grounding wire with an appropriate grounding plug. The plug must be plugged into an outlet that is properly installed and grounded in accordance with all local codes and ordinances. If power code is not equipped (option) then appropriate grounding shall be provided upon installation.

 **Warning**

Improper installation of the grounding plug is able to result in a risk of electric shock. When repair or replacement of the cord or plug is required, do not connect the grounding wire to either flat blade terminal. The wire with insulation having an outer surface that is green with or without yellow stripes is the grounding wire.

 **Note**

Check with a qualified electrician or serviceman when the grounding instructions are not completely understood, or when in doubt as to whether the product is properly grounded. Do not modify the plug provided; if it does not fit the outlet, have the proper outlet installed by a qualified electrician.

 **Note**

This pump must be connected to a grounded, metallic, permanent wiring system, or an equipment-grounding terminal or lead on the product.

Warning

Do not connect an input power source to a thermal sign takeoff wire. A motor is damaged.

3.5 Fluctuations in the power voltage and frequency

Standard: Rotation electricity machine general rules

JIS C 4034-1:1999, JEC-2137-2000

To the voltage change and frequency change in Domain A, in main rated values, it operates continuously, and can be used practically convenient, and to the voltage change and frequency change in Domain B, it shall operate with main rated values and shall be used practically convenient.

However, operation with "it is convenient and safe is maintained on "practical use, it means not resulting in the grade which shortens a life remarkably, and the characteristic, a temperature rise, etc. do not apply correspondingly in the state of rating. Moreover, main rating shows rated torque (N·m).

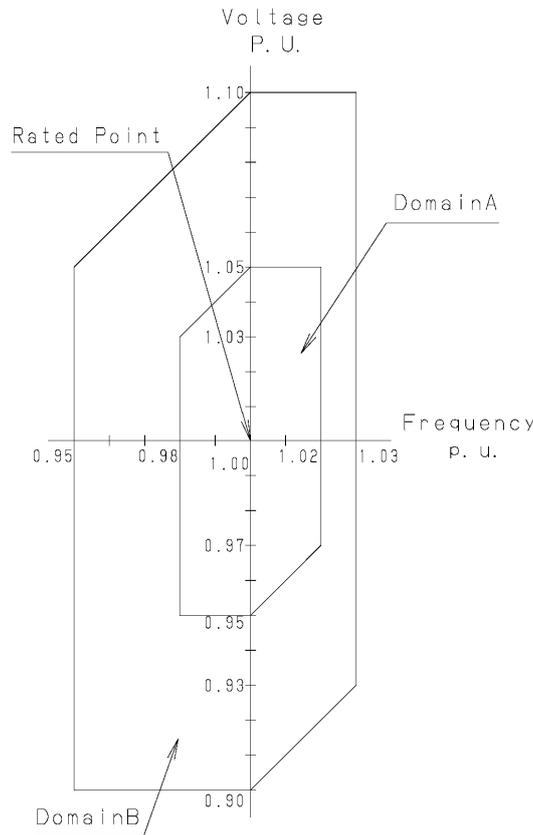


Fig. 7 Change region of the voltage and frequency

 **Warning**

Before connecting wires, turn off the power switch. Never perform wiring with the power supplied as an electric shock will occur. Connect the earth wire correctly. Failure to do so may result in electric shock if a failure or earth leakage occurs. Installation of a dedicated earth leakage breaker is also recommended.

 **Warning**

A start mode of this motor is condenser start / operating method. The switching of the start / operation reaches by electronic governor switch. When I apply voltage of the inverter control because I use an electron element for electronic governor switch, malfunction of the electronic governor switch occurs and causes the motor stumbling-block. Please never apply voltage of the inverter control.

 **Caution**

Perform electric wiring correctly in accordance with the “Electric Equipment Technical Standard” and “Internal Wiring Regulation.” Incorrect wiring will result in fire.

 **Caution**

Install an overload protector suitable for the capacity of the motor. If an overload protector is not installed, or if an overload protector that is unsuitable for the motor capacity is installed, the motor will be damaged leading to fire.

4. Operation

4.1 Cautions for operation

Warning

There is a risk of explosion. Never block the outlet or operate the pump with equipment mounted at the outlet side which blocks the passage of gas. Otherwise, the pump internal pressure increases causing the pump to explode, the oil level gauge to protrude or the motor to be overloaded.

This pump is not resistant to pressure. The internal pump pressure is limited to 0.03 MPa (gauge pressure).

Caution

Do not stop and restart this equipment repeatedly. When restarting this equipment, make sure that the pump (motor) has stopped completely before turning on the power. If the pump (motor) has not stopped completely and the equipment is restarted, the pump's (motor) current rises and either the protection device is activated or the motor may become damaged.

Note

- ① In the process of manufacturing semiconductors, pump oil may deteriorate over a very short period of time. It is recommended that the pump oil should be replaced within 10 days after starting use of the pump, and the replacement frequency of the pump oil should be decided based on the contamination level of the pump oil.
- ② If the pump evacuate a lot of moisture, replace the oil frequently. If the pump is used with gas which contains a lot of moisture, the lubricity of the pump oil deteriorates and corrosion of the pump's components advance, causing the pump to malfunction.
- ③ If chemicals including acid has been evacuated, the pump may become rusty while it is not being operated (i.e. overnight), making operation impossible. If such chemicals are evacuated, replace the pump oil immediately.
- ④ Solvents which deteriorate the lubricity of the pump oil will cause scoring, etc. If such a solvent is evacuated, replace the oil.
- ⑤ If operation is performed continuously at a high evacuation pressure of 1 kPa or more, a large amount of pump oil is consumed, causing a shortage of oil and insufficient lubrication of the pump. If such a condition continues, components will rapidly wear and become scored. Avoid continuous operation at a high evacuation pressure as much as possible and, without fail, add pump oil.
- ⑥ This pump uses carbon in internal member of framework. Oil may become black when it runs. However, there is not it in the aberration.

4.2 Start of operation

To start operation, close leak valve (B), open vacuum valve (A) to the inlet port, and turn on the power switch. Then the pump starts beings to exhaust (see Fig. 5).

Caution

- ① Firing hazard. Do not touch the motor or the pump body as these surface temperature should be higher (70~80°C).
- ② If operation is performed at high pressure, oil mist is generated at the outlet side. Install an oil mist trap or connect a duct to discharge the oil mist outside the room. Or, install a ventilator.

Note

When the pump does not rotate correctly, take the following measures.

- a) Check the amount of oil, and adjust if necessary.
- b) In an environment where the ambient temperature is low, if the pump is left unused for a long time (three days or longer), the pump oil enters the cylinder. (This phenomena cannot be avoided even if the pump pressure is released to atmospheric pressure after last using the pump.) If the pump is restarted in this condition, an overload is applied to the pump and the overload protector may actuate. In such a case, turn the pump on and off several times in short intervals.

Note

After operating the pump for several hours, oil temperature inside the pump may be risen up to 70~80°C. If the temperature is over this, the pump may have troubles. Maintain the pump, or contact us.

4.3 Stopping the operation

To stop operation, close vacuum valve (A), open leak valve (B) quickly, and turn the power switch off (see Fig. 5).

Caution

Do not stop and restart this equipment repeatedly. When restarting this equipment, make sure that the pump (motor) has stopped completely before turning on the power. If the pump (motor) has not stopped completely and the equipment is restarted, the pump's (motor) current rises and either the protection device is activated or the motor may become damaged.

Caution

Firing hazard. As the surface of the pump body is higher temperature (70~80°C) during operation, do not touch it until cooled down after stopping the pump.

4.4 Backflow preventer (G-101S)

A backflow preventer is incorporated into the pump to prevent the oil from flowing back while the pump is stopped.

The backflow preventer actuates in the case of an emergency including power failure. So, after the pump is stopped due to a power failure, follow the procedures mentioned in “4.3 Stopping the operation” to stop the operation.

Note

- ① To stop the pump, always close vacuum valve (A) and then open leak valve (B). If this procedure is neglected, the pump oil fills the cylinder, making restart difficult or causing damage to the pump. The pump oil also may flow back to the vacuum chamber side.
- ② If vacuum valve (A) is not closed, air may leak into the device side through the pump increasing the pressure.

4.5 Thermal protector

A manually reset thermal protector is incorporated in the motor in order to interrupt the power circuit of the motor and prevent damage to the motor when an over current flows through the motor due to a stop in rotation or overload resulting from the pump malfunctioning during operation. The type of thermal protector to be used depends on the pump model.

Table 3 Characteristics of the thermal protector

Model		Operating characteristic	
		Operation temperature	The temperature that can reset
G-101D, G-101S	1 ϕ , 100V	100°C \pm 7°C	74°C \pm 7°C
	1 ϕ , 200V	100°C \pm 7°C	74°C \pm 7°C

When thermal protector operates, at first I disconnect the plug, and please inform our company. The motor is very hot when the thermal protector has actuated. Never touch it with your hand. When the cause of the malfunction has been eliminated, check that the motor has cooled down, and then press the manual reset button to restart operation (see “6.5 Trouble check list”).

Caution

Firing hazard. As the surface of the pump body is higher temperature (70~80°C) during operation, do not touch it until cooled down after stopping the pump.

4.6 Gas ballast valve (G-101D)

The pump is equipped with a gas ballast valve in order to evacuate vapor and condensable gases such as solvent vapor.

Evacuated condensable gas that liquefies in the compression and pressurization processes of the pump is mixed with the pump oil and starts circulating through the pump together with the oil. In such a case, the same effect as when oil of a high steam pressure is used is produced, and the ultimate pressure of the pump increases. Moreover, the lubricity of oil deteriorates and the service life of the shaft seal is shortened.

If air or dry nitrogen enters through the gas ballast valve just before the compression and

pressurization processes of the pump, condensable gas will not liquefy and will be exhausted together with air through the outlet valve. When the gas ballast valve is used, the “gas ballast effect” increases as the pump temperature becomes high. So, before evacuating condensable gas, perform operation for approximately 20 minutes with the gas ballast open, and after the pump temperature reaches approximately 50 ~ 65°C, open vacuum valve (A) and continue operation. If the temperature is low, a satisfactory “gas ballast effect” is not achieved.

If the gas ballast valve is left open when condensable gas is not evacuated, not only does the pump oil scatter and power is lost, but also the ultimate pressure increases. Furthermore, since the gas ballast valve’s capacity to process condensable gas is limited, condensable gas remains in the pump oil when a lot of condensable gas is exhausted or when condensable gas (air and gas containing small amounts of moisture and other vapor which make the oil dirty) is exhausted without opening the gas ballast valve. In such a case, perform non-load operation with vacuum valve (A) closed and the gas ballast valve open. Then the oil temperature increases and the pump oil is purified due to the effect of the gas ballast valve. Continue non-load operation with the gas ballast valve closed until the specified pressure is reached. If the pump oil is not cleaned even a long time, replace the pump oil.



Caution

As the surface of the pump body is high temperature (70~80°C) during operation, do not touch it besides the gas-ballast valve.
Start operation after closing the gas-ballast valve.



Note

If the gas ballast valve is left open without condensable gas being exhausted, the pump oil scatters, power is lost, or the ultimate pressure increases.
Close the gas ballast valve when condensable gas is not exhausted.

4.7 Operation in cold climates

4.7.1 Operational method

In winter-season, when operating the pump in cold environment , it will be difficult to start up the pump. It is called over-load operation because of high viscosity of the pump oil at low temperature.

If the pump does not restart, warm up the pump oil, or try to keep the pump within the operating temperature for a little while and then turn on the power again.

When the pump stops after rotating for a few seconds, open leak valve (B) and continuous operation may become possible. After the pump has warmed up, close leak valve (B) and return to ordinary operation.

4.7.2 Important points

This pump uses the motor of the condenser start / operation type.

Start / the operation performs switching in electronic governor switch made guts in motor.

When start of the pump is difficult, pump does a restart action for approximately five seconds, when it cannot start, thermal protector operates for conservation.

When thermal protector operated, I disconnect the plug, and let me reset thermal protector (see “4.5 thermal protector ”).

4.8 An inlet port, exhaust port replacement (an option)

4.8.1 The switching method of the inlet port

When I change an inlet port of the pump, please choose 40□ flange of our optional inlet port. The switching method excludes screw fixing an inlet port, and please attach an inlet port to change.

4.8.2 The switching method of the exhaust port

When I change exhaust port of the pump, please choose G1 size screw of our optional exhaust port. The switching method removes exhaust port fixing, and please attach an exhaust port to change.

4.9 Installation of the oil mist trap (Option)

An oil mist trap can be installed in order to remove oil mist from the pump. As such an oil mist trap, models OMT-100A and OMI-100 are available. Remove the standard exhaust port from the exhaust port of the pump and install an oil mist trap instead. The oil mist trap not only prevents oil mist generation but also reduces exhaust noise.

For details, refer to the instruction manual for the oil mist trap.

4.10 Restriction on operation when the oil mist trap is installed

When using the oil mist trap, there are some restrictions on operation. When the filter is clogged, replace it.

The internal pump pressure is limited to 0.03 MPa (gauge pressure). When the pressure measured at the outlet side has reached 0.03 MPa (gauge pressure), replace the oil mist trap filter.



Warning

Be sure to observe the restrictions on operation when the oil mist trap is installed. There is a risk of explosion. When the filter is clogged, replace it.

5. Pump Performance

5.1 Ultimate pressure

The term “ultimate pressure” as employed in the catalogue and in this manual is defined as “the minimum pressure obtained by the pump without the introduction of gas from the pump inlet (i.e. the non-load condition).” For this pump, measurement is performed using the specified pump oil with only a Pirani vacuum gauge installed at the pump inlet port.

Note that the Pirani gauge shows values approximately five to ten times higher than those shown by the McLeod gauge. This is because condensable gas components (mainly moisture) included in the measured air are removed when the McLeod gauge is used.

Also, the actual ultimate pressure of the vacuum device becomes higher than that noted in the catalogue for the following reasons.

- (1) The vacuum gauge is installed at a distance from the pump, and the steam and a variety of gases are generated by water droplets and rust on the inside walls of the pump and piping.
- (2) Gasifying of volatile components which have dissolved in the pump oil. (Deterioration of pump oil)
- (3) Existence of a gas supply source including vacuum leakage in the vacuum path.

5.2 Pumping speed

The pumping speed of the pump depends on the type and pressure of the gas to be evacuated. The pumping speed usually reaches the maximum at a high pressure range, and it gradually decreases as the pressure reduces.

The nominal pumping speed of this pump is the maximum pumping speed when dry air is evacuated. Fig. 6 shows the relationship between the evacuation pressure and pumping speed.

5.3 Power requirement

The power required to operate the pump is the total of the power required to overcome the rotational resistance of the pump (mechanical work) and the power required to compress the air (compression work), and reaches a maximum at an inlet evacuation pressure of around 2.7×10^4 to 4×10^4 Pa. If the inlet evacuation pressure has reduced to 13.3 Pa or less, the compression work is considerably reduced and more power is consumed in mechanical work.

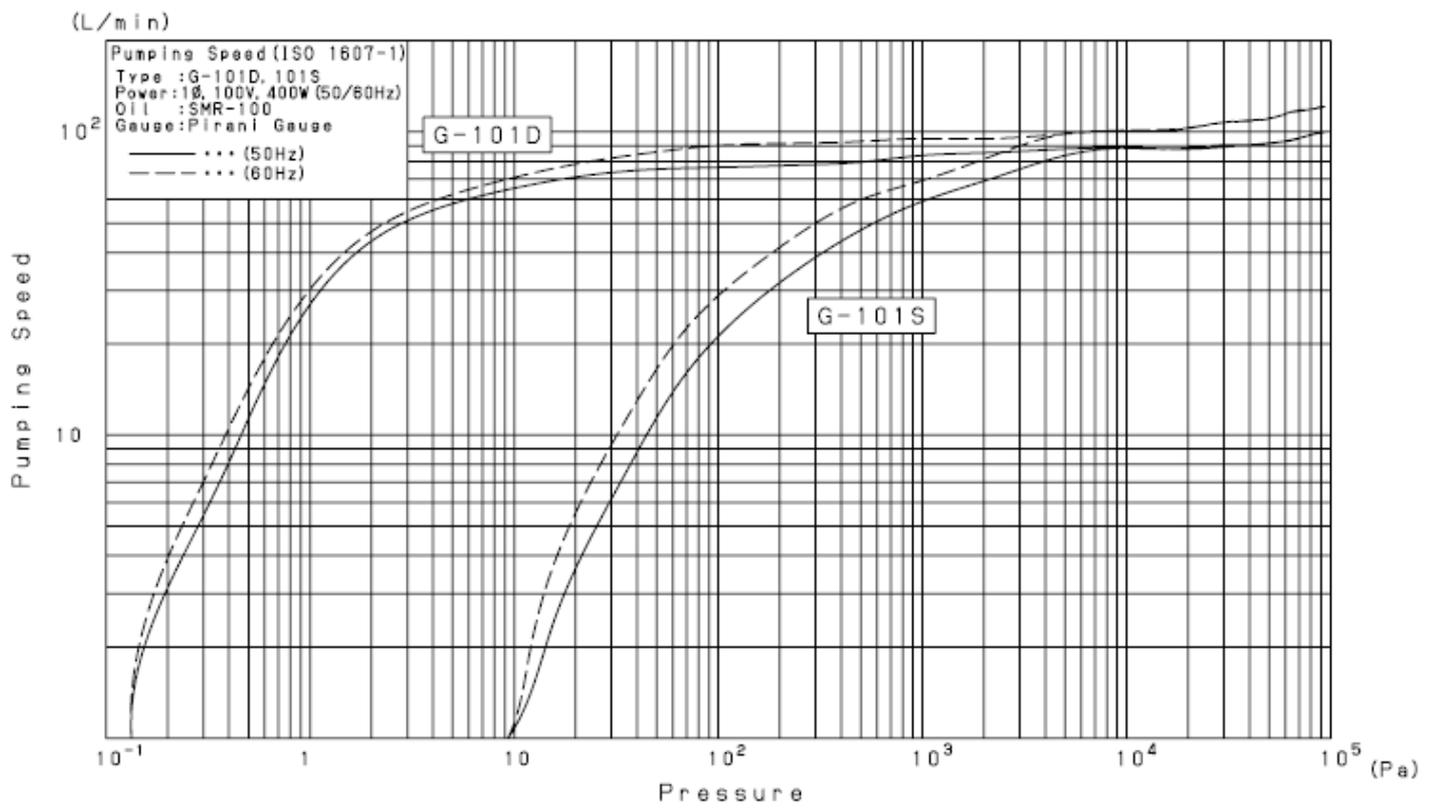


Fig. 8 Pumping speed curve

6. Maintenance, Inspection and Repair

6.1 Maintenance

Check the following during operation at least once every three days.

- (1) Amount of pump oil (To be within the range shown with red lines on the oil level gauge)
- (2) Discoloration of the pump oil
- (3) Abnormal sound
- (4) Problem with the motor current value
- (5) Oil leak from the oil seal

If there is any problem, take proper measures in accordance with “6.5 Trouble check list.”

6.2 Periodic inspection

The items to be checked should be changed as necessary depending on the environment where the pump is used. However, always check the following in order to prevent a malfunction and to lengthen the service life of the pump.



Caution

- ① Turn off the power before starting inspection and do not turn it on while inspection is in progress. Doing so will result in injury.
- ② The pump is very hot immediately after it is stopped. Wait for a while until the pump has cooled down completely and then start inspection. There is a risk of burns.

1) Periodic replacement of the pump oil

The pump oil deteriorates with operation. Check the viscosity and level of contamination of the pump oil with the oil level gauge, and replace the pump oil in good time. If the pump oil is replaced periodically, the deterioration of the pump oil is minimized and the service life of the pump is lengthened.

If operation is continued with a lot of moisture mixed with the pump oil, the ultimate pressure will not reach the standard value, the movement at the section where the mechanical friction is generated becomes slow, and the pump finally becomes damaged. Replace the pump oil in accordance with “6.3 Replacement of the pump oil.”

Table 4 Periodic inspection table

Frequency	Item	Details	Measures
Once/3 days	Oil	Amount	Refill the oil.
		Color (Reddish brown, dark blown, and cloudy white are not good.) This pump uses carbon in internal member of framework. Oil may become black when it runs. However, there is not it in the aberration.	Replace the oil.
	Sound	Abnormal sound	Check nuts and bolts for looseness. If not clear, contact us.
	Vibration	Abnormal vibration	
	Current value	Difference from the rated value	Check the cause of an overload. If not clear, contact us.
Once/week	Surface temperature	Surface temperature (The temperature higher than the room temperature by 45°C or more is abnormal.)	Check the cause of an overload. If not clear, contact us.
	Oil leakage	Oil leakage from the shaft seal section and plugs.	Replace seals, or contact us.
Once/3,000 operation hours or once/6 months	Inlet filter	Clogged with dust	Clean the inlet filter.
	Oil	Even if no problem is recognized, be sure to replace the oil.	Replace the oil.
Once/year	Coupling	Damage or fracture	Replace the coupling.
	Oil pump filter	Clogged with dust	Clean the oil pump filter.

2) Inspection of the amount of pump oil

Refill the pump oil so that the pump oil level is kept within the range of the red lines showing the upper and lower limits on the oil level gauge during operation.

3) Inspection of oil leakage

When oil leaks from the shaft seal section or drain plug seal section, repair is required. Our specified O-rings and seals are always available from the service departments shown at the back of this manual. When necessary, contact them.

4) Inspection of inlet filter

If the inlet filter is clogged with dust included in the evacuated gas, the pump's efficiency may deteriorate.

5) Inspection of abnormal sounds and vibration

Check the nuts and bolts for looseness.

6) Inspection of the coupling

Check the coupling which connects the main pump unit and motor of the pump for damage. If cracks or fractures are found on the coupling, replace it in accordance with “6.4 Replacement of the coupling.”

7) Inspection of the oil mist trap

When using the oil mist trap in replacement of the standard exhaust port, pay attention to the clogging of the filter in the oil mist trap. If the clogging advances, evacuated gas cannot be exhausted any longer, which causes the oil gauge to protrude and oil leakage from the shaft seal section or drain plug seal section. The maximum internal pump pressure is 0.03 MPa (gauge pressure).

When the pump is operated continuously for a long time or when the pump is extremely contaminated with evacuated gas, overhaul is required. Contact the nearest sales or service department among those listed at the back of this manual.

 **Danger**

When requesting the manufacturer's service department to overhaul the pump, always write the type of the vacuumed gas on the “Pump Usage Check Sheet” attached at the back of this manual and submit it. Note that if toxic gases are exhausted, both the pump itself and pump oil will become contaminated. Please be sufficiently aware that use with some gases will preclude overhaul.

6.3 Replacement of the pump oil

The pressure of the vacuum device may increase due to the deterioration of the pump oil. In such a case, close the inlet port of the pump and check that the specified ultimate pressure has been reached. If not, replace the pump oil. If substances having a high vapor pressure (such as moisture or solvents) are mixed with the pump oil, or if sludge is accumulated at the bottom of the pump, the ultimate pressure cannot be reached with only one replacement and the pump oil must be replaced several times. The deterioration of the pump oil is caused not only by the contamination due to evacuated gas but also by the changes in the properties of the pump oil itself (depending on the operation time). Periodic replacement in accordance with Table 4 showing an oil replacement guide is recommended.

 **Danger**

Keep in mind that if the pump is used for exhausting toxic gas, both the main pump unit and pump oil will become contaminated.

 **Caution**

- ① Wear protective equipment such as rubber gloves and safety goggles.
- ② Be sure to read the attached “Safety Data Sheet” before adding oil. If the oil accidentally comes into contact with your hands or enters your eyes, take proper measures in accordance with the section “First-aid treatment” shown in “Safety Data Sheet”

 **Note**

Use only oils specified by us. If other oils are used, the pump performance will deteriorate or its life will shorten.

< Pump oil replacement procedure >

- (1) Release the pump inlet port to the atmosphere and operate the pump for five seconds. The oil remaining in the pump is discharged efficiently.
- (2) Remove the exhaust port and drain plug to discharge the pump oil.
- (3) Mount the drain plug, and add the required amount of the new specified pump oil through the lubrication port (see Fig. 4).
- (4) If the pump oil is contaminated extremely, add new pump oil and perform operation for a while (several minutes) to clean the pump. Repeat this a few times.
- (5) After replacing with the new pump oil, operate the pump and when the pump has become warm, check the ultimate pressure.
- (6) If the pump oil is so dirty that oil sludge accumulates at the bottom of the pump, the specified ultimate pressure even after the pump oil is replaced. In such a case, overhaul the pump.

6.4 Replacement of the coupling

A coupling is used at the section connecting the pump main unit and the motor. It is recommended that this coupling be periodically inspected once a year or so. If the corner is chipped or cracked, replace it. If the pump is started and stopped hundreds of times a day, increase the inspection frequency.

To take out the coupling, discharge the pump oil, and remove the hex. socket set bolt (which fixes the coupling to the motor shaft from the back of the base) and the four bolts (which fix the pump main unit to the base). Then the coupling can be taken out to remove the bolt which fixes the coupling by the side of casing. After inspecting the coupling, re-attachment should attach coupling so that the plane part of a casing side rotor shaft and a stop screw with the hex. socket set bolt, and please perform it in the way shown in Fig. 9.

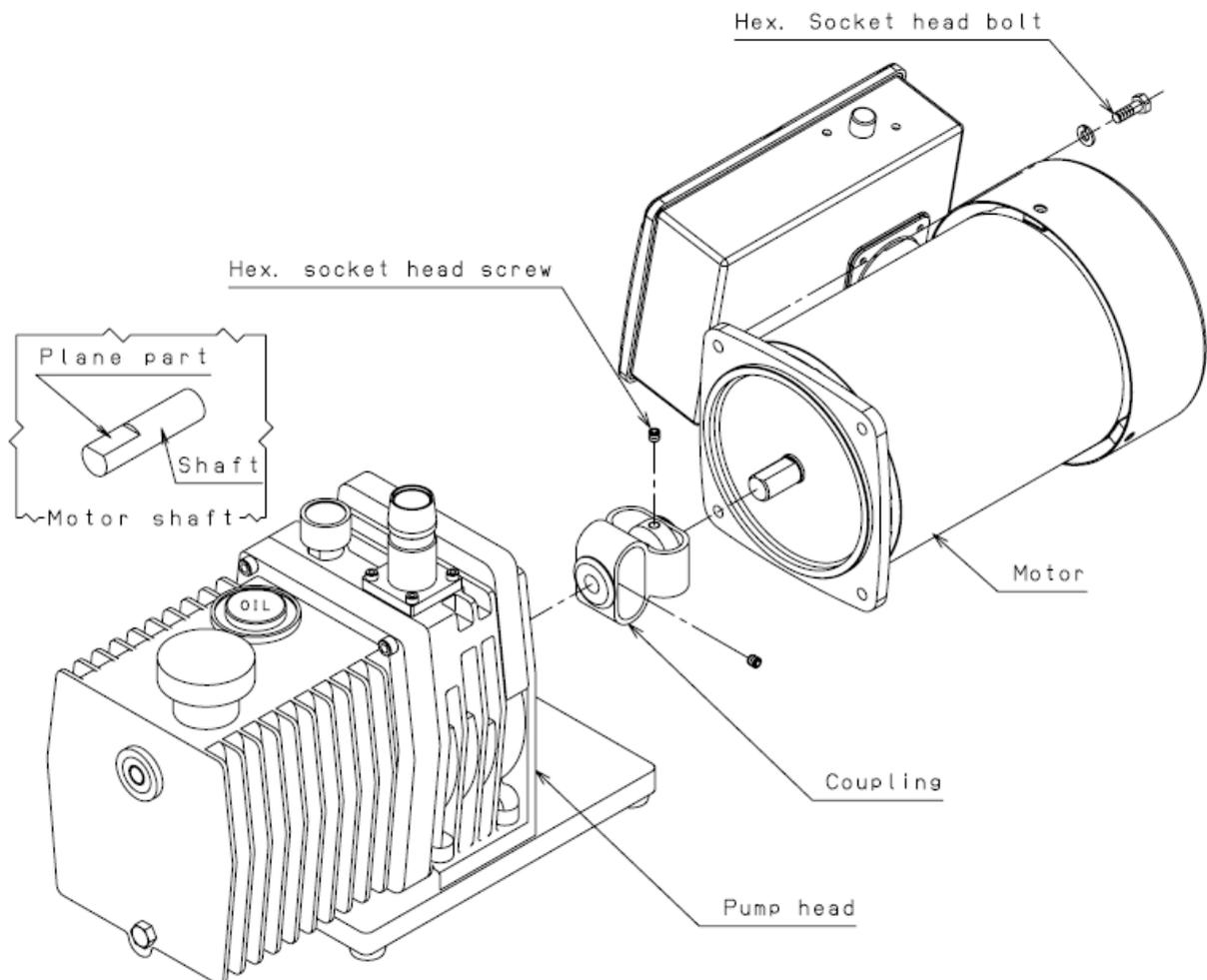


Fig. 9 Replacement of the coupling

Connect the concave section (female) of the pump unit with the convex section (male) of the motor, push the motor into the pump so that both connecting surfaces come completely into contact with each other, and fix the motor with bolts.

6.5 Trouble check list

Table 5 Trouble check list

Problem	Cause	Measures	Reference
The pump does not rotate.	① The pump is not connected to the power supply.	① Connect the pump to the power supply.	3.4
	② The power switch is not turned on.	② Turn on the power switch.	4.2
	③ Problem with power supply voltage	③ Set the power supply voltage to within $\pm 10\%$ of the rated voltage.	3.5
	④ The overload protector has actuated.	④ Disconnect the plug and push the reset button. Press the reset button.	4.6
	⑤ The motor malfunctions.	⑤ Replace the motor.	
	⑥ Low ambient temperature has increased the oil viscosity.	⑥ Increase the ambient temperature to 7°C or more.	4.4
	⑦ The entrance of foreign matter into the pump caused the rotor to burn out.	⑦ Overhaul (replace the cylinder and rotor).	6.2
	⑧ Moisture or solvents were sucked in, forming rust inside the pump.	⑧ Overhaul (replace the cylinder and rotor).	6.2
	⑨ Reaction product accumulated in the pump when the pump stops after exhausting reactive gas.	⑨ Overhaul (clean the pump inside and remove reaction products).	
	⑩ Components inside the pump have burnt out.	⑩ Overhaul (replace the damaged components).	
The pump's rotation is unstable.	① Problem with power supply voltage	① Set the power supply voltage to within $\pm 10\%$ of the rated voltage.	3.5
	② Defective wiring to the pump	② Perform wiring to the pump again.	3.4
	③ Low ambient temperature has increased the oil viscosity.	③ Increase the ambient temperature to 7°C or more.	4.4
	④ Foreign matter has entered the pump.	④ Disassemble and clean the pump to eliminate foreign matter.	
The pressure does not decrease.	① The pump is too small for the volume of the vacuum chamber.	① Select another pump.	5.2
	② The pressure measurement method is not correct.	② Measure the pressure correctly.	5.1
	③ The vacuum gauge is not suitable.	③ Measure with a calibrated vacuum gauge suitable for the pressure range.	
	④ The pipe connected to the inlet port is small, or the piping distance is long.	④ Use pipes having a diameter larger than the inlet port diameter, or reduce the distance from the vacuum chamber.	5.1
	⑤ The wire mesh at the inlet port is clogged.	⑤ Remove the piping from the upper section of the inlet port, and clean the wire mesh.	6.2

Problem	Cause	Measures	Reference
The pressure does not decrease.	⑥ The specified amount of oil has not been added.	⑥ Add the specified amount of oil.	3.2
	⑦ The oil has deteriorated.	⑦ Replace the oil.	6.3
	⑧ Leakage occurs from the pipe connected to the pump.	⑧ Locate the leakage with a leakage detector and stop the leakage.	6.3
	⑨ Our specified oil is not being used.	⑨ Overhaul the pump and replace with oil specified by us	
	⑩ Oil does not circulate, or the oil hole of the cover is clogged.	⑩ Overhaul and clean the oil hole.	
Abnormal sound is generated.	① Problem with power supply voltage	① Set the power supply voltage to within $\pm 10\%$ of the rated voltage.	3.5
	② The motor malfunctions.	② Replace the motor.	3.2
	③ Foreign matter has entered the pump.	③ Eliminate the foreign matter and overhaul the pump.	
	④ The specified amount of oil has not been added.	④ Add the specified amount of oil.	
	⑤ Oil does not circulate, or the oil hole of the cover is clogged.	⑤ Overhaul and clean the oil hole.	
	⑥ Components inside the pump have burnt out.	⑥ Overhaul (replace the damaged components).	
Pump surfaces are extremely hot (50 °C or more higher than the room temperature)	① Continuous operation at high evacuation pressure	① A pump is suspended or suction pressure is set to 1kPa or less.	3.2
	② The specified amount of oil has not been added. (If the oil amount is not sufficient, the cooling effect of the pump will be reduced.)	② Add the specified amount of oil.	
	③ The temperature of the evacuated gas is high.	③ Mount cooling equipment such as a gas cooler at the inlet side.	
	④ Oil does not circulate, or the oil hole of the cover is clogged.	④ Overhaul and clean the oil hole.	
A lot of oil splashes out from the exhaust port.	① The pump is been filled in excess of the specified amount.	① Discharge the oil until it reduces to the specified amount.	3.2
	② Continuous operation is performed at a high evacuation pressure.	② Install an oil mist trap at the outlet side.	4.8
The oil leaks outside the pump.	① Deterioration of the O-ring and the oil seal of the case and cover	① Check and replace the O-ring and oil seal.	6.2

7. Disposal

Follow state law and local government regulations for disposal of the pump.



Caution

- ① When a harmful toxic gas has been exhausted, ask a specialist for waste disposal. Not only the pump itself but also the pump oil become toxic.
- ② For the disposal of pump oil, follow the instructions given under “Cautions for disposal” in “Safety Data Sheet.”

8. Maintenance parts

8.1 Maintenance parts list

Table 6 G-101D Maintenance parts list

Product name	No.	Parts name	O'ty
G-101D Maintenance kit A	17	Outlet valve	4
	18	Outlet valve spring	4
	38	O-ring_S-6	1
	39	O-ring_S-24	1
	40	O-ring_S-55	1
	41	O-ring_P-8	1
	42	O-ring_P-12	1
	43	O-ring_P-24	1
	44	O-ring_V-160	1
	45	Oil seal_HTC-17-40-9	1
	46	Oil seal_SCY-17-30-7	1
	47	Oil seal_SC-15-30-7	1
	70	O-ring_P-20	1
G-101D Maintenance kit B	6	1st vane	2
	7	Vane spring	8
	12	2nd vane	2
	14	Middle pin	4
	28	Oil level gauge	1
	31	Coupling $\phi 14 \times \phi 16$	1
	36	Bearing_6203	1
	37	Bearing_6002	1
	-	Tube	1
	-	Sleeve_SN1/8	2

Note : For the relationship between components, see the disassembly drawing.

Table 7 G-101S Maintenance parts list

Product name	No.	Parts name	O'ty
G-101S Maintenance kit A	11	Outlet valve	3
	12	Outlet valve spring	3
	41	O-ring_S-4	1
	42	O-ring_S-6	1
	43	O-ring_S-18	2
	44	O-ring_S-24	1
	45	O-ring_S-55	1
	46	O-ring_P-8	1
	48	O-ring_P-24	1
	49	O-ring_V-160	1
	50	Oil seal_HTC-17-40-9	1
	51	Oil seal_SCY-17-30-7	1
	52	Oil seal_VC-15-30-5	1
	79	O-ring_P-20	1
G-101S Maintenance kit B	6	Vane	2
	7	Vane spring	4
	8	Middle pin	2
	31	Oil level gauge	1
	34	Coupling $\phi 14 \times \phi 16$	1
	39	Bearing_6203	1
	40	Bearing_6002	1

Note : For the relationship between components, see the disassembly drawing.

8.2 Disassembly drawing

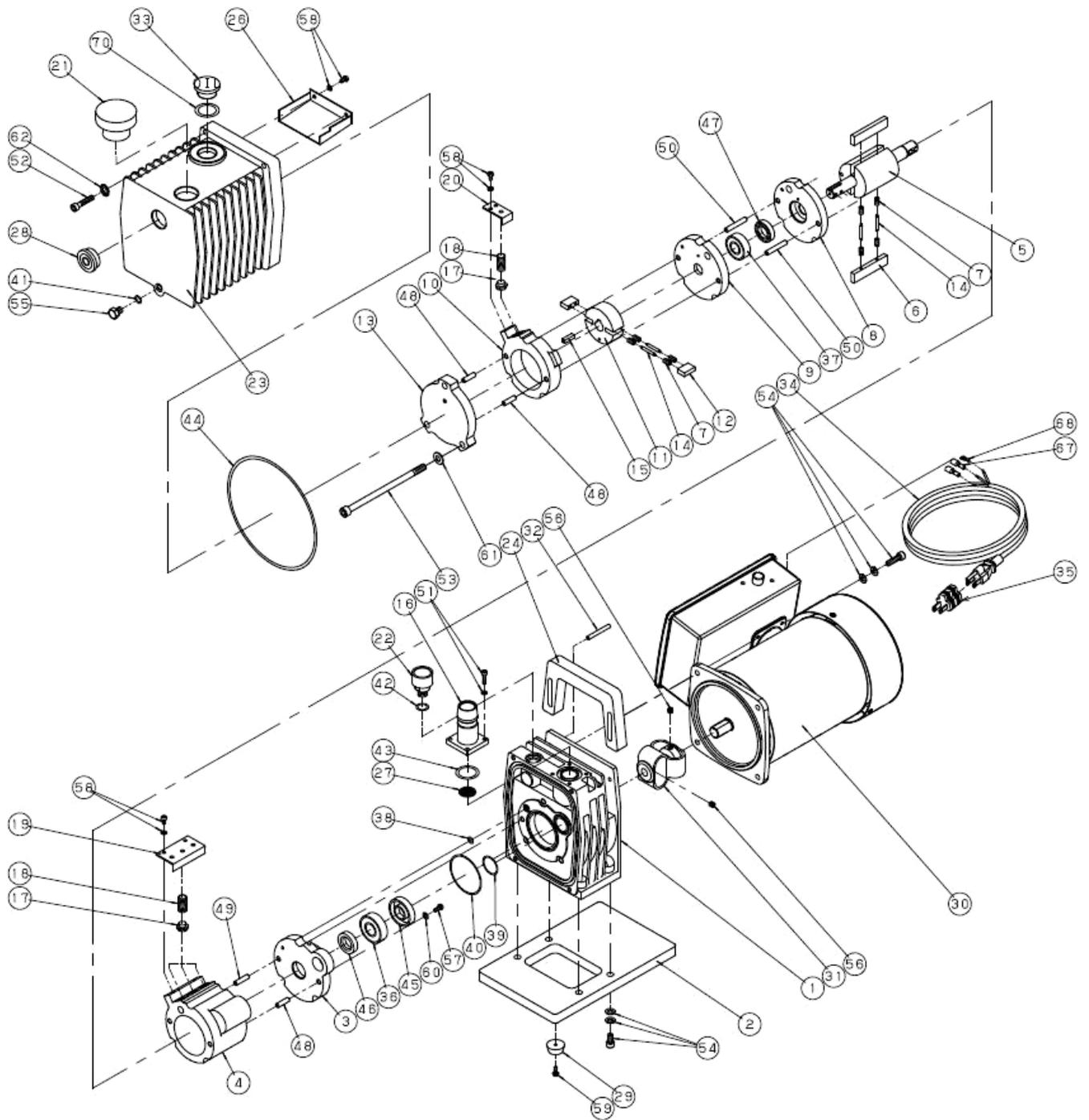


Fig. 10 Disassembly drawing of G-101D oil sealed rotary vacuum pump

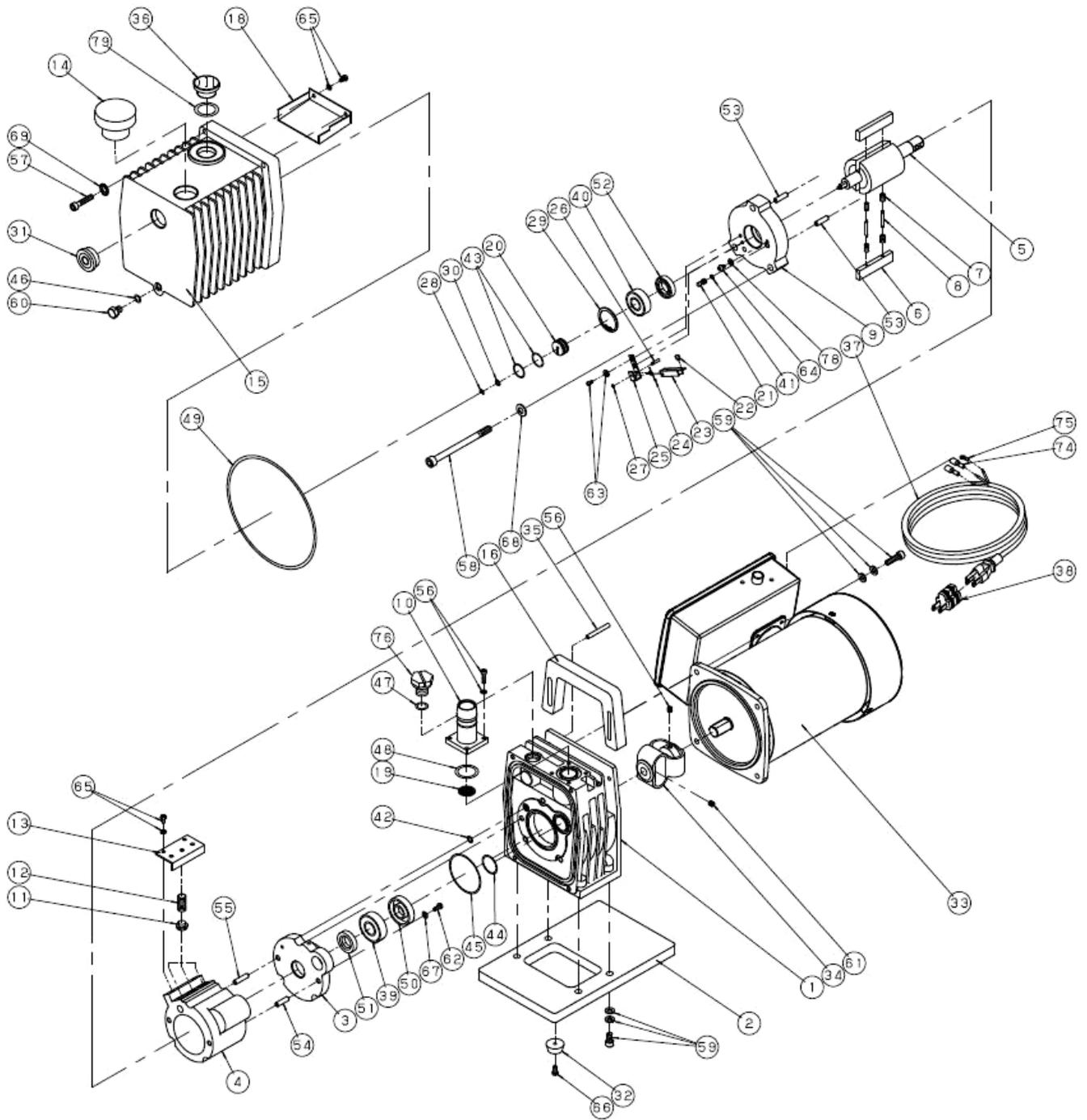


Fig. 11 Disassembly drawing of G-101S oil sealed rotary vacuum pump

Warranty

- (1) The warranty for this pump (this equipment) extends for a period of one year from the date of shipment.
- (2) Any malfunctions or defects which occur under normal usage conditions during the warranty period will be repaired free of charge.

Note, the warranty stated here is an individual warranty covering the pump. In addition, the scope of the warranty coverage concerning repairs is limited to the repair and/or replacement of parts.

Normal usage conditions refer to the following:

- a) Ambient temperature and humidity during operation: 7 - 40°C, below 85% RH
- b) Type and temperature of the exhausted gas: Dry air or dry nitrogen, 7 ~ 40□
- c) Operation in accordance with the user manual

- (3) Repair fees will incur during the warranty period for the following cases:

- a) Malfunctions due to a natural disaster or fire.
- b) Malfunctions caused by special atmospheric conditions, such as salt damage, inflammable gas, corrosive gas, radiation or pollution.
- c) Malfunctions caused by usage conditions that differ from those stated in the user manual (performance specifications, maintenance and inspection, etc.).
- d) Malfunctions caused by modifications or repairs carried out by a party other than the manufacturer, or by a service company not approved by the manufacturer.
- e) Malfunctions caused by noise (electric disturbance).
- f) Malfunctions that occur when not using a rated power supply.
- g) Malfunctions that occur when there is an abnormal rise in internal pressure due to the pump exhaust outlet being blocked during operation, etc.
- h) Malfunctions that occur, when the pump is damaged as a result of being dropped or falling, etc.
- i) Malfunctions which are determined by the manufacturer's technical personnel to be caused by conditions that do not comply with the usage conditions for this vacuum pump.
- j) Malfunctions due to the replacement of consumables.

- (4) Disclaimer

- a) We shall not be liable for any malfunctions of our products caused by the customer, regardless if the malfunction does not fall within the warranty period, nor shall we be liable for any loss of opportunity for the customer's clients or for compensation for any damages to other products, labor costs, production loss, transportation expenses and other related work.
- b) We shall not be liable for any claims and patent infringements, including secondary damages, filed a claim by a third party against the customer.

Usage Status Check Sheet (for use in Instruction Manual)

* For the purpose of safety control of repair personnel, fill in within the heavy line frame and attach the sheet to the item of which repair is requested.

* In case this sheet were not attached or filled in, your request of repair and service may not be accepted.

* In accordance with the Private Information Protection Law, the provided information will be used only for determining the cause of failure and whether detoxifying washing should be conducted. It will never be provided to any third person.

Model Name: _____ Manufacturer's Serial No.: _____

1. Inhaled Gas * Please be sure to fill in.

(1) Whether there is harmful effect on human bodies Yes No (Sing your name below.)

(2) Whether there is unusual smell Yes No

(3) Type and Name of Gas: _____

* Industrial Safety and Health Law designates particular substances as the materials to be notified.

2. Usage Status

Operation Method: Approx. () hours per day, () years and () months

Continuous Operation Intermittent Operation

Usage: _____

3. Failure Status Unusual Noise Abnormal Pressure Abnormal Actuation

Oil Leakage Other Symptoms: _____

4. Detail of Request Repair (Overhaul) Regular Checks

5. Others: _____

Company Name: _____ Personnel in charge: _____

Address: _____

Tel: _____ Fax: _____ E-mail: _____

Agent Name: _____ Personnel in charge: _____

Address: _____

Tel: _____ Fax: _____

* In case you do not have any direct transaction with us, please be sure to fill in the agent name.

6. Confirmation

The gas and substance used in this pump or unit is harmless to human bodies, or it is not contaminated by any substance harmful to human bodies.

Signed _____ (seal) Date: _____

* In order to avoid a trouble during transportation, please evacuate oil from any oil pump before shipping.

* You are requested to ship the package to our Service Division (CS Center). (See the attached list of addresses.)