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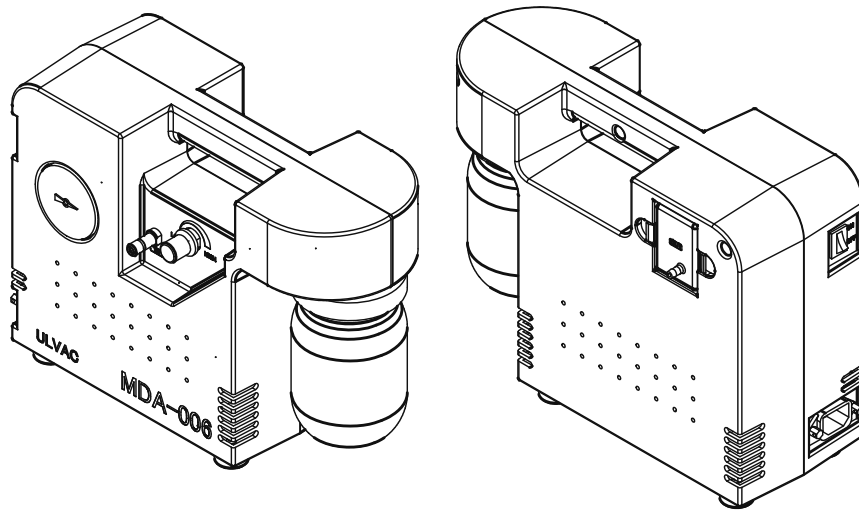
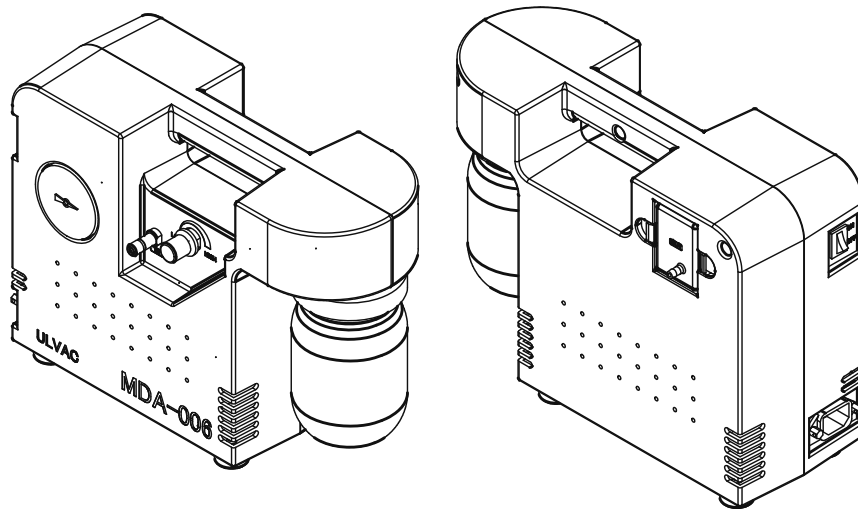
Instruction Manual  
For  
Portable Aspirator

Model Name

MDA-006

**Request to Users**

Please read this manual thoroughly to ensure safe and effective use of the equipment.

Keep this manual in a place where it can be referred to at any time and look it carefully.

The contents of this instruction manual are subject to change without prior notice due to improvements in performance and the functions of the product

# Table of contents

Items with shade include description on safety

(0) To safely use this equipment	0 1
1. Power supply	0 2
2. Environment	0 3
3. Setting	0 3
4. Operation	0 4
5. About maintenance, repair, the disposal	0 5
(1) Before using	1
1. Target users	1
2. Read the manual thoroughly	1
3. This manual in a safe place	1
4. Statutory requirements for disposal	1
5. Safety during repair	1
(2) Product overview	2
1. Purpose of this product and actions that are prohibited	2
2. Safety device and its purpose and functions	2
3. Product specification	3
4. Individual unit specifications	3
5. Standard accessories	4
6. Switch and valve usage methods	4
7. Switch and valve layout	5
(3) Opening the package and installation	6
1. General cautions	6
2. Package upon delivery	6
3. Installation site	6
4. Power supply	7
5. Power supply cord fixation metal fittings	8
(4) Operations	9
1. Risks and safety measures upon operations	9
2. Portable aspirator operation procedure	9
2-1 Preparation	9
2-2 Operations	9

2-3	Stop	1 0
3.	Measures upon abnormal conditions	1 0
3-1	During power outages	1 0
3-2	Operation after recovery from power outage	1 0
(5)	Maintenance and repair	1 1
1.	Risks and safety measure upon maintenance and repair	1 1
2.	Maintenance and repair that can be made by user	1 1
3.	Removal, maintenance, and installation of devices	1 2
3-1	Exhaust tank	1 2
4.	Maintenance and inspection points	1 3
5.	Troubleshooting	1 3
6.	Storage of equipment	1 4
7.	Consumable list	1 4
(6)	Disposal	1 4
1.	Cautions upon disposal	1 4
(7)	Conclusion	1 4

Warranty

Use situation check seat

Business, the service sector and the address

## (O) To safely use this equipment

Thank you for purchasing this product. This instruction manual is a guidebook to have used this device safely. The basic instructions and control point necessary for the handling and the check maintenance point are listed. It is well use after a reading by the contents which explain in this instruction manual to prevent a serious accident. As for the instruction manual and this safe guide, we hold copyright or other. Intellectual property rights. Please do not reproduce some or all instruction manuals without our permission. And a correction or the delivery cannot do it.

Read this section before using the MDA-006.

Follow the instructions below to safely use the device and prevent personal injuries from occurring.

Please comply with them all the time.

The meanings of the figure sign of the text are as follows.

 Danger	Incorrect handling of the equipment is very likely to result in death or serious injury to the operator.
 Warning	Incorrect handling of the equipment may result in death or serious injury to the operator.
 Caution	Incorrect handling of the equipment may result in light or medium injuries to the operator or damage to the equipment.
	This indicates action or practice that should be made.
	Always make connection with the earth
	This indicates the action or practice that should be prohibited.
	Do not disassemble.
	Do not touch.

# 1. Power supply

 <p>Warning</p>	 <p>Voltage confirmation</p>	<p>Please do not use it other than the product rating voltage. A thermal protector does not operate normally, It causes the damage by a fire of the motor, the fire.</p>
	 <p>Check the cable capacity</p>	<p>Avoid using an extension cable. However, if it must be used by necessity, use the following cable. For 100-V power supply: 1.25 mm<sup>2</sup> or more If a thinner cable is used, this may cause overheating, ignition, or fire.</p>
	 <p>Avoid this action</p>	<p>Do not perform the desorption of the power supply cord by a wet hand It causes the electric shock.</p>
	 <p>Avoid this action</p>	<p>Do not use with the primary side cable bundled up. In addition, do not place any objects on the cable. Otherwise, such action may cause electrical shock or fire.</p>
	 <p>Check the ground</p>	<p>Connect the primary side cable to a grounded outlet. If using a plug adapter, securely connect to ground. There is a risk of electrical shock when damaged or if there is a short circuit.</p>

## 2. Environment

 <b>Danger</b>	 Avoid this action	Absorption gas leaks from other than the inlet port of this device, a vent and the inside. Please do not absorb a poisonous gas.
	 Avoid this action	This product does not have explosion-proof design, and thus use in environments where inflammable substances are present should be avoided. Otherwise, explosion could occur, causing fire and burns.
 <b>Warning</b>	 Avoid this action	The pump is extremely hot during operation. Do not place flammable objects around it. This can cause fire.
	 Environmental confirmation	Please do not put a thing disturbing the ventilation around a device. There might be a burn due to the abnormal fever, the fire.
 <b>Caution</b>	 Environmental confirmation	Pump is heated during operation. Please do not put a thing affected by heat in the neighborhood.

## 3. Setting

 <b>Warning</b>	 Environmental confirmation	Install the equipment where the following conditions are satisfied. <ul style="list-style-type: none"> <li>a) Flat surface</li> <li>b) Floor with sufficient strength</li> <li>c) Well-ventilated place</li> <li>d) Place without direct sunlight</li> <li>e) Room with temperatures between 7° C and 40° C.</li> <li>f) Location where there is no risk of fire</li> <li>g) Place without electrical noises, which may cause adverse effect to the product.</li> <li>h) Without oil mist and the grease spot.</li> <li>i) Without dew condensation and dust.</li> </ul> Otherwise, operation failure or durability degradation may occur.
---	---	--

## 4. Operation

 Danger	 Open exhaust outlet	<p>Make sure the exhaust outlet is open.</p> <p>Please remove a thing disturbing the passage of the gas of the vent side.</p> <p>The maximum internal pressure is 0.03 MPa (gauge pressure).</p>
	 Avoid this action	<p>This product is not a medical treatment apparatus. Because it causes an injury and the accident, please do not use it for the human body.</p>
 Caution	 Avoid this action	<p>Please do not breathe in a material unsuitable for EPDM/FPM rubber. (including a liquid, the gas including corrosive gas, an organic solvent, a liquid and condensability gas, mineral oil, steam, the sugar)</p> <p>When absorb it, it causes the decline of the driving abnormal cause and durability.</p>
	 Avoid this action	<p>Please do not absorb the solid body.</p> <p>The solid is clogged up by the internal plumbing. (Symptom :) The peculiar noise / flow quantity is short ;).</p>

## 5. About maintenance, repair, the disposal

 <b>Warning</b>	 Resolution action	<p>Please do not perform the resolution, the disassembly of a resin cover and the exhaust tank. It causes the trouble.</p>
 <b>Caution</b>	 Maintenance and repair range	<p>Please confirm the following item before driving.</p> <ol style="list-style-type: none"> <li>1) Is nothing in the collection jar?</li> <li>2) Does not an abnormal sound occur?</li> <li>3) Do not block up the vent?</li> </ol>
	 Wear protective equipment	<p>Wear protective equipment (gloves, glasses, etc.) when draining fluid from inside the collection jar or replacing the diaphragm. Some aspirated gases may cause toxic substances to be left in the collected liquid and inside the pump.</p>
	 Comply with regulations	<p>The device and all accessory parts must be disposed of in accordance with legal and regional municipality regulations. Comply with regulations upon disposal. For details on disposal, contact us.</p>
	 Avoid this action	<p>Do not use other manufacturer's optional parts for modification. We do not assume any responsibility for any damage due to such modifications.</p>

## ( 1 ) Before using

### 1. Target users

Only persons who have used aspirators or been trained based on this manual may operate this equipment.

### 2. Read the manual thoroughly

Please read this manual thoroughly in order to use the equipment in a safe and correct manner.

Please pay particular attention when reading the section “To Safely Use This Equipment” .

### 3. Keep this manual in a safe place

Keep this manual in a safe place.

After reading this manual, be sure to keep it in a safe place where it is readily accessible to other users.

### 4. Statutory requirements for disposal

Dispose of the equipment parts, please obey of all laws and ordinances and a local government rule.

Please obey a rule for disposal.

### 5. Safety during repair

When requesting repairs to this product, provide a full description of the conditions of use (particularly any use of dangerous materials) for the safety of repair personnel.

In this case, fill in the Use Condition Check Sheet and attach to the product.

If the use conditions are unknown, repair may be refused.

## (2) Product overview

### 1. Purpose of this product and actions that are prohibited

This device is a portable aspirator equipped with a diaphragm type vacuum pump. This is different from a water aspirator in that the ultimate pressure it can reach is not affected by tap water pressure or temperature and can be used even in locations with no tap water source.

Adjust it by any pressure and can use it while watching the vacuum gauge which put on a panel.

To safely and correctly use this product, avoid the following actions that are prohibited.

 Avoid this action	<ul style="list-style-type: none"> <li>▪ This device is only for vacuum exhaust. Please do not run for long time in the atmospheric pressure neighborhood. It causes the trouble.</li> <li>▪ This product is not a medical treatment apparatus. Because it causes an injury and the accident, please do not use it for the human body.</li> <li>▪ Please do not use this device for pressurization use. (Please refrain from the absorption of the air which pressurized.)</li> <li>▪ Please do not do the resale of this device, the resolution, remodeling.</li> </ul>
--	--

### 2. Safety device and its purpose and functions

Item	Purpose	Functions
 Avoid this action	Diaphragm type dry vacuum pump protection	The diaphragm type dry vacuum pump circuit protector operates to stop power flow.
Operation with the safety devices above disabled is prohibited.		

### 3. Product specifications

Item	Specifications
Pressure adjustment range	6.6 ~ 98 kPa (absolute pressure)
Pressure indication	Analog (Bourdon tube vacuum gauge)
Aspiration pressure adjustment	Needle valve (VAC. CONT)
Aspiration port external diameter	φ 9 mm
Vent outer diameter	φ 6 mm
Collection jar	450 cc × 1
Vacuum pump	DA-15D (Special)
Structure	Diaphragm type dry vacuum pump
Pumping speed	6 L/min (50 Hz)
Ultimate pressure	6.6 kPa (absolute pressure)
Ambient pressure for usage	7 ~ 40 °C
Power Supply	AC 100 V 0.5 A or above (50 Hz) 0.5 A or above (60 Hz)
Power requirement	0.05 kVA
Dimensions	137.5 mm(W) × 313 mm(D) × 241 mm(H)
Weight	Approx. 4 kg

### 4. Individual unit specifications

Unit	Model and specifications	Quantity
1) Diaphragm type dry vacuum pump	<ul style="list-style-type: none"> <li>• Model DAP-6D (Special)</li> <li>• Pumping speed 6 L/min (50Hz)</li> <li>• Ultimate pressure 6.6 kPa (absolute pressure)</li> <li>• Required power supply 0.05 kVA</li> </ul>	1 set
2) Bourdon tube vacuum gauge	<ul style="list-style-type: none"> <li>• Model TE-22SP-10mm</li> <li>• Pressure graduation 0-0.1 MPa (absolute pressure)</li> </ul>	1 pc.

## 5. Standard accessories

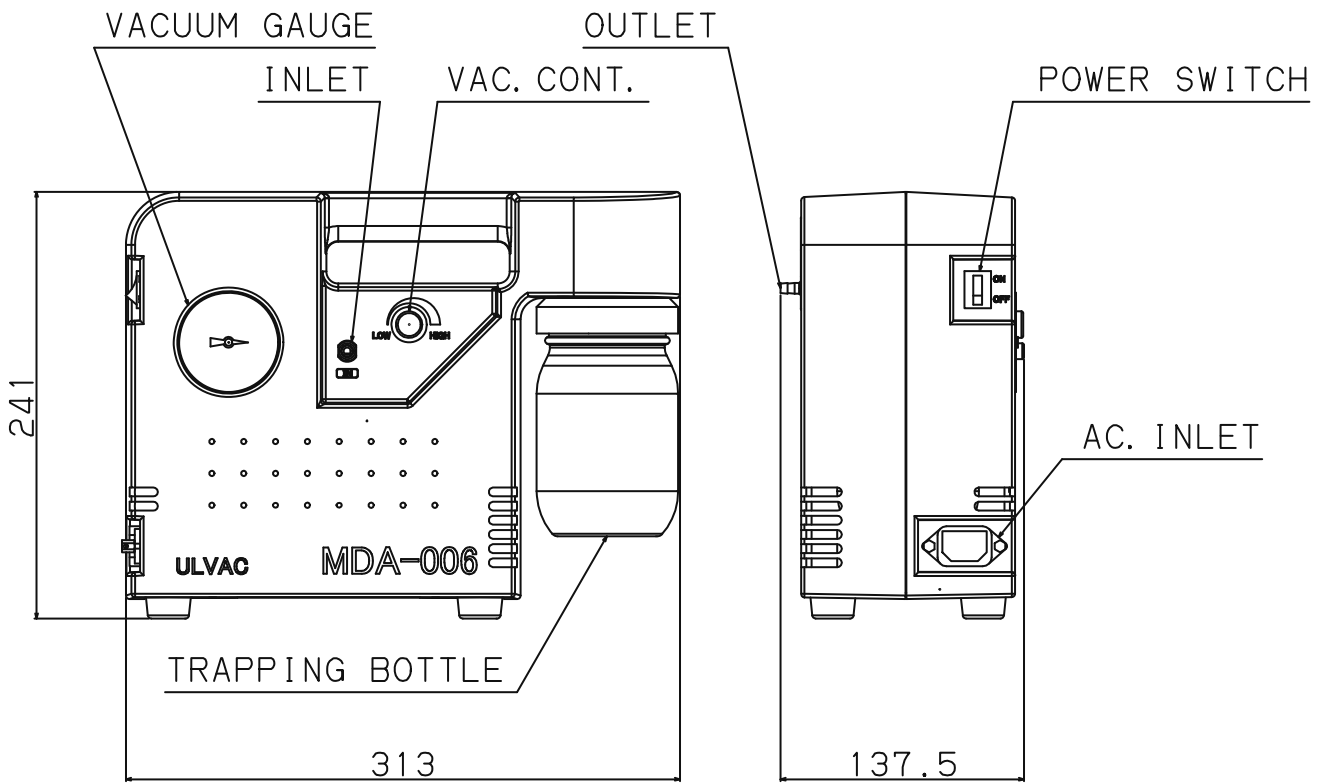
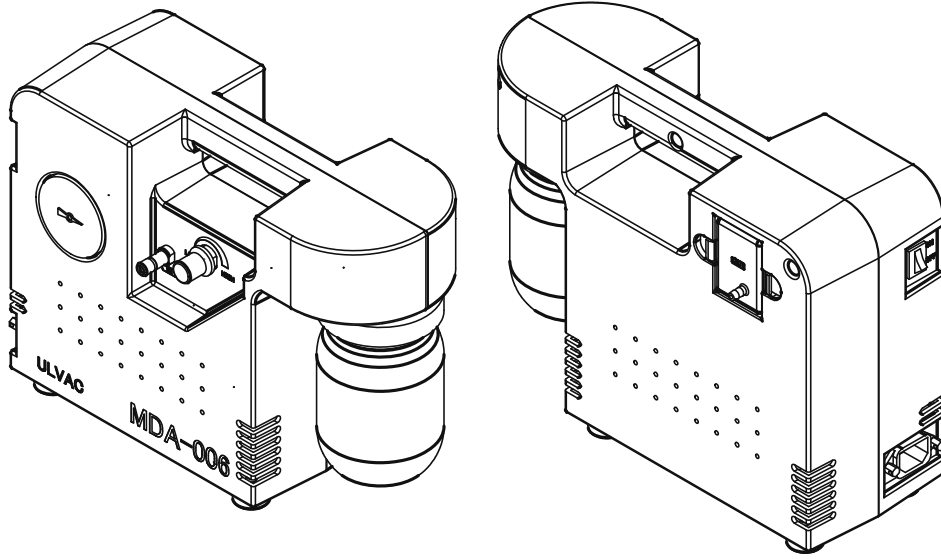
1) Power supply cable	- 100 V single phase, with power outlet, 2 m	1 pc.
2) Power supply cord fixation metal appliance	- 100V power supply cord use	1 pc.
3) Plug adapter	- Plug with grounding wire for 100-V power supply	1 pc.
4) Aspiration port cap	- for $\phi 9$	1 pc.
5) Instruction Manual	- This book	1 pc.

## 6. Switch and valve usage methods

 Check	<p>The switch becoming OFF before an operation start, Valve (VAC.CONT) please start operation in confirming that LOW (shut) becomes in a state.</p>
--	---

Name	How to Operate
Power switch	Snap switch Manual ON-OFF
Needle valve (V A C C O N T . )	Screw tightening type HIGH: Clockwise (open) LOW: Counter clockwise (close)

## 7. Switch and valve layout (VAC. CONT)

### (3) Opening the package and installation

#### 1. General cautions

 <p>Check</p>	<ol style="list-style-type: none"><li>1) Is the product what you requested?</li><li>2) Are the accessories and all necessary parts included?</li><li>3) Check that no damage has occurred during shipping.</li><li>4) Maintain a space of more than 15 cm from the equipment for safety upon installation.</li></ol>
--	--

#### 2. Package upon delivery

The main unit is shipped packed in a cardboard box.

#### 3. Installation site

 <p>Warning</p>	 <p>Check the environment</p>	<p>Install the equipment where the following conditions are satisfied.</p> <ol style="list-style-type: none"><li>a) Flat surface</li><li>b) Floor with sufficient strength</li><li>c) Well-ventilated place</li><li>d) Place without direct sunlight</li><li>e) Room with temperatures between 7° C and 40° C.</li><li>f) Location where there is no risk of fire</li><li>g) Place without electrical noises, which may cause adverse effect to the product.</li><li>h) Without oil mist and the grease spot.</li><li>i) Without dew condensation and dust.</li></ol> <p>Otherwise, operation failure or durability degradation may occur.</p>
 <p>Check</p>	<p>After the installation, confirm that the product has clearance of at least 15cm from the wall for safety</p>	

## 4. Power supply

 Warning	 Check the capacity	Please prepare the following: Single-phase 100 V (50Hz/60Hz) A thermal protector does not operate other than the rating voltage normally and causes the damage by a fire of the motor, the fire.
Primary side power supply scramble specifications (part attached to the device) Power supply cord 2,000mm in length for 100V 3P plug (with an adapter) with tip ground		
 Warning	 Check the ground	Connect the primary side cable to a grounded outlet. If using a plug adapter, securely connect to ground. There is a risk of electrical shock when damaged or if there is a short circuit.
	 Check the cable capacity	Avoid using an extension cable. However, if it must be used by necessity, use the following cable. For 100-V power supply: 1.25 mm <sup>2</sup> or more If a thinner cable is used, this may cause overheating, ignition, or fire.
	 Avoid electrical shock	When connect power supply cord, connected to the device earlier, and please insert it in a power supply after having fixed it by power supply cord fixation metal appliance. It causes the electric shock.
	 Avoid this action	Do not use with the primary side cable bundled up. In addition, do not place any objects on the cable. Otherwise, such action may cause electrical shock or fire.

## 5. Power supply cord fixation metal fittings

Warning

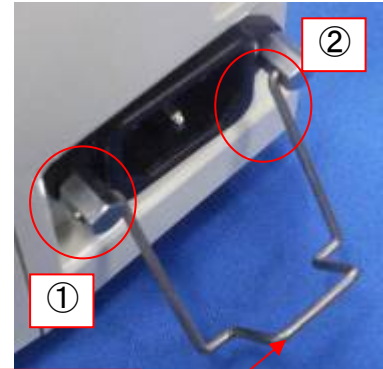
Check

Please fix power supply cord by power supply cord fixation metal fittings by all means.

### Power supply cord fixation metal fittings installation procedure

Please attach the power supply cord fixation metal fittings according to the following procedures.

- ① Please hang one of power supply cord fixation metal fittings on the hole of the power supply cord plug mouth (inlet) side from the inside.
- ② Please hang it on the hole of the power supply cord plug mouth (inlet) side from the inside while pushing the other side of power supply cord fixation metal fittings.


Power supply cord fixation metal appliance

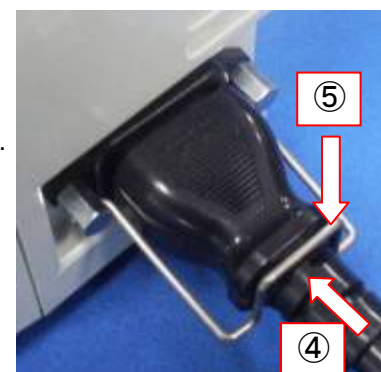- ③ Lift power supply cord fixation metal fittings.

- ④ Insert power supply cord well.
- ⑤ Lower the power supply cord fixation metal fittings and fix power supply cord.

With the above installation of power supply cord fixation metal fittings is completion.

## (4) Operations

### 1. Risks and safety measures upon operations

 Danger	 Open exhaust outlet	Make sure the exhaust outlet is open. Please remove a thing disturbing the passage of the gas of the vent side. The maximum internal pressure is 0.03 MPa (gauge pressure).
 Caution	 Avoid this action	Please do not breathe in a material unsuitable for EPDM/FPM rubber. (including a liquid, the gas including corrosive gas, an organic solvent, a liquid and condensability gas, mineral oil, steam, the sugar) When absorb it, it causes the decline of the driving abnormal cause and durability.

### 2. Portable aspirator operation procedure

<p>2-1 Preparation</p> <ol style="list-style-type: none"><li>1) Open a packing case and take the internal cover, and please take out a device having Toride of the device.</li><li>2) Please put a device in the installation location.</li><li>3) Please the outskirts of the device keep away a device from the wall at least 15cm to cool it.</li><li>4) A cap is attached to an inlet port. Remove a cap, and please connect a pipe (including the rubber pipe). (please make the vent opening.)</li><li>5) Set the VAC CONT. valve to LOW.</li><li>6) Please insert power supply cord in the main body of the device. Please attach a power supply cord fixation metal appliance. Please check the attached power supply cord fixation metal appliance. Please insert a power supply cord plug in the nearest outlet. And please confirm that do not have the slack.</li></ol> <p>This concludes operation preparation.</p>
<p>2-2 Operations</p> <ol style="list-style-type: none"><li>1) Please turn on POWER SWITCH. Start the diaphragm vacuum pump.</li><li>2) Set the VAC CONT. to HIGH. Adjust pressure to desired pressure. (Setting the pressure to 60 kPa or above may cause the Bourdon tube vacuum gauge needle to sway slightly, however this is not abnormal. The swaying is caused by the diaphragm pulsing.)</li></ol>

### 2 – 3 Stop

- 1) Turn the POWER OFF.  
Stop the diaphragm vacuum pump.
- 2) Turn knob (VAC. CONT) for a direction of HIGH, and please cancel a vacuum.
- 3) Remove a liquid in the disassembly with a collection jar (Turn it clockwise and remove it)

### Precautions

- 1) If stopped while there is still a vacuum, the pump will not start when the device is restarted.  
Turn the VAC. CONT knob to the right to let air in a remove the vacuum.
- 2) If liquids collect to approximately half of the collection jar, stop a device and, with a rubber cap, turn a slow bottle counterclockwise, and please remove it. And please throw away a liquid.
- 3) This product has an automatic return type thermal protector built-in. When motor temperature rises by a stop and overload by the trouble of the product while driving, cut off the power supply circuit of the motor automatically, and this prevents the damage by a fire of the motor.  
When a thermal protector operates, at first turn off a switch, and please contact us. The motor neighborhood (main body center) please never touches it by hand then because temperature rises.  
If it is removed a cause of the trouble, please reopen driving after having confirmed that main body temperature falls.
- 4) Aspirating mists will have little effect. (Some types of mist may decrease performance.) Directly aspirating liquids will cause the pump to lock and trip the circuit protector.  
Aspirating mists will have little effect. (Some types of mist may decrease performance.) Directly aspirating liquids will cause the pump to lock and trip the circuit protector.
- 5) Please do not breathe in a material unsuitable for EPDM/FPM rubber. (including a liquid, the gas including corrosive gas, an organic solvent, a liquid and condensability gas, mineral oil, steam, the sugar)  
When absorb it, it causes the decline of the driving abnormal cause and durability.

## 3. Measures upon abnormal conditions

### 3 – 1 During power outages

- 1) Turn the POWER OFF,
- 2) Set the VAC CONT. to HIGH.

### 3 – 2 Operation after recovery from power outage

Follow the steps "2-1 Preparation" and "2-2 Operations" under 2. Portable Aspirator Operation Procedure".

## (5) Maintenance and repair

### 1. Risks and safety measure upon maintenance and repair

 <p>Caution</p>	 <p>Wear protective equipment</p>	<p>Wear protective equipment (gloves, glasses, etc.) when draining fluid from inside the collection jar or replacing the diaphragm.</p> <p>Some aspirated gases may cause toxic substances to be left in the collected liquid and inside the pump.</p>
--	--	--

### 2. Maintenance and repair that can be made by user

<p>1) Disassembly of the collection jar</p> <p>2) Cleaning of the collection jar</p> <p>Please do not perform the resolution except the above, the repair</p>
---

### 3. Removal, maintenance, and installation of devices

 <b>Caution</b>	 Wear protective equipment	Please wear a tool for protection (gloves, glasses) for the cleaning of the exhaust tank. Some aspirated gases may cause toxic substances to be left in the collected liquid and inside the pump.
	 Avoid this action	Please do not remove the cover of the exhaust tank, and the main body of tank and the cover are glued together in a liquid gasket. A liquid gasket comes off when take off a cover, and a leak occurs than the tank which attached again.

3-1 The liquid disposal is the following methods installation, disassembly Please go.

1) Disassembly order

- Put a finger in a knob and push it inward.

- Pull it slowly toward.

- Because a silicon tube is connected to the tank back, please outrun. (please be careful so that a tube is not in the main body inside then)

2) Liquid disposal

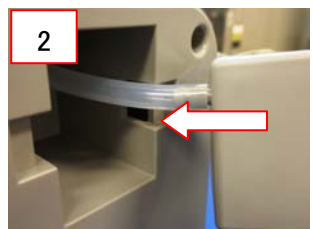
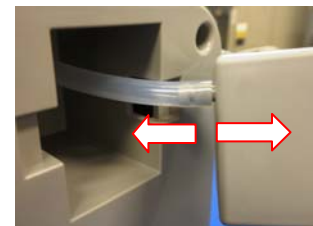
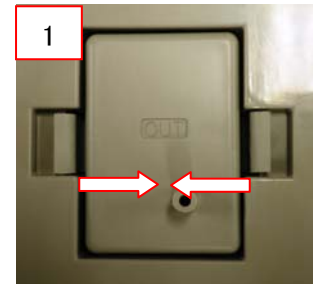
- Do not take the cover of the tank and Slant it and discard a liquid. Do not make reverse and verticality. For the lower half of the cover, and the liquid gasket applies

3) Installation order

- Connect a tube to a tank.

- Push a tank. Please push the tank slowly. A tube might be destroyed.

- The part of the nail please push it until a sound sounds with a click.

#### 4. Maintenance and inspection points

Unit	Maintenance and Inspection	Timing of Maintenance and Inspection
Aspirator	Verify there is no liquid in the collection jar.	Before every use.
	Dispose of liquid in the collection jar every time the device is stopped.	After every use.
	Please throw away the liquid in the exhaust tank.	After every use.
	Verify performance (ultimate pressure).	As necessary

#### 5. Troubleshooting

Symptom	Cause	Troubleshooting
Does not start Stop that is running	The operation of the thermal protector due to the abnormal fever	Switch it off, and please remove a trouble cause
	Power supply abnormality	Please confirm whether an appropriate power supply is supplied
	Vacuum start	Please perform atmosphere opening
Ultimate pressure is poor, ultimate pressure is unstable or Evacuation performance is low.	Alien substance mixture to the plumbing	Please remove an alien substance Alien substance mixture please do not use it on a possible condition
	Diaphragm type vacuum pump is defective.	Refer to the "Diaphragm Type Vacuum Pump User's Manual".
	The ambient temperature at the installation site is high.	Please use it in the range of atmosphere temperature 7 degrees Celsius - 40 degrees Celsius
	Valve operation is not appropriate	Perform opening and shutting confirmation of valve (VAC. CONT), and please adjust it to appropriate pressure
	The slack of the collection jar	Please attach a sampling bottle appropriately
	Water was in the pump	Switch on this device and adjust pressure to half atmospheric pressure, and please confirm the pressure after driving for approximately 30 minutes
	Failure in measuring instrument	Replace the measuring instrument.

When a symptom is not improved even if give the measures mentioned above, please contact us

## 6. Storage of Equipment

Observe the following guidelines to store the equipment.
<p>1) Place of storage</p> <ul style="list-style-type: none"> <li>- Floor with sufficient strength</li> <li>- Location with good ventilation</li> <li>- Location with no vibration.</li> <li>- Location without direct sunlight</li> <li>- Location where room temperature is below normal temperature.</li> <li>- Location where no chemicals or gases or other substances which can have a corrosive effect on the exterior of the device are present.</li> </ul>
<p>2) Cautions before and upon storing</p> <ul style="list-style-type: none"> <li>- Attach a cap to an inlet pipe.</li> <li>- A collection jar makes an open state.</li> <li>- Discard the liquid in the exhaust tank.</li> </ul>

## 7. Consumable List

Location of Use	Parts	Specifications	Material	Quantity
Rack	Rubber foot	----	NBR	4
	Collection jar	450cc	Soda-lime glass	1

※ When seem to be trouble, please ask the dealer for repair after communication without disintegrating in own.

### (6) Disposal

#### 1. Cautions upon disposal

 <b>Caution</b>	 Comply with regulations	<p>The device and all accessory parts must be disposed of in accordance with legal and regional municipality regulations.</p> <p>Comply with regulations upon disposal.</p> <p>For details on disposal, contact us.</p>
Applied regulations: Rules of waste disposal and cleaning		
Disposal methods: 1) Transportation: Industrial waste contractors 2) Disposal: Industrial waste contractors		

### (7) Conclusion

<p>Explained the handling, the driving method of this aspirator, but the content is limited to a general matter as things mentioned above. If there are a questionable point and a trouble, ask to have or we talk about communication.</p>
---

## Warranty

- (1) The warranty for this pump (this **aspirator**) extends for a period of one year from the date of shipment.
- (2) Any malfunctions or defects which occur under normal usage conditions during the warranty period will be repaired free of charge.

Note, the warranty stated here is an individual warranty covering the pump. In addition, the scope of the warranty coverage concerning repairs is limited to the repair and/or replacement of parts.

Normal usage conditions refer to the following:

- a) Ambient temperature and humidity during operation: 7 - 40°C, below 85% RH
- b) Operation in accordance with the user manual

- (3) Repair fees will incur during the warranty period for the following cases:

- a) Malfunctions due to a natural disaster or fire.
- b) Malfunctions caused by special atmospheric conditions, such as salt damage, inflammable gas, corrosive gas, radiation or pollution.
- c) Malfunctions caused by usage conditions that differ from those stated in the user manual (performance specifications, maintenance and inspection, etc.).
- d) Malfunctions caused by modifications or repairs carried out by a party other than the manufacturer, or by a service company not approved by the manufacturer.
- e) Malfunctions caused by noise (electric disturbance).
- f) Malfunctions that occur when not using a rated power supply.
- g) Malfunctions that occur when there is an abnormal rise in internal pressure due to the pump exhaust outlet being blocked during operation, etc.
- h) Malfunctions that occur, when the pump is damaged as a result of being dropped or falling, etc.
- i) Malfunctions which are determined by the manufacturer's technical personnel to be caused by conditions that do not comply with the usage conditions for this vacuum pump.
- j) Malfunctions due to the replacement of consumables.

- (4) Disclaimer

- a) We shall not be liable for any malfunctions of our products caused by the customer, regardless if the malfunction does not fall within the warranty period, nor shall we be liable for any loss of opportunity for the customer's clients or for compensation for any damages to other products, labor costs, production loss, transportation expenses and other related work.
- b) We shall not be liable for any claims and patent infringements, including secondary damages, filed a claim by a third party against the customer.

Usage Status Check Sheet (for use in Operation Manual)

- \* For the purpose of safety control of repair personnel, fill in within the heavy line frame and attach the sheet to the item of which repair is requested.
- \* In case this sheet were not attached or filled in, your request of repair and service may not be accepted.
- \* In accordance with the Private Information Protection Law, the provided information will be used only for determining the cause of failure and whether detoxifying washing should be conducted. It will never be provided to any third person.

Model Name: \_\_\_\_\_ Manufacturer's Serial No.: \_\_\_\_\_

1. Inhaled Gas \* Please be sure to fill in.

(1) Whether there is harmful effect on human bodies Yes No (Sing your name below.)

(2) Whether there is unusual smell Yes No

(3) Type and Name of Gas: \_\_\_\_\_

\* Industrial Safety and Health Law designates particular substances as the materials to be notified.

2. Usage Status

Operation Method: Approx. ( ) hours per day, ( ) years and ( ) months

Continuous Operation  Intermittent Operation

Usage: \_\_\_\_\_

3. Failure Status  Unusual Noise  Abnormal Pressure  Abnormal Actuation

Oil Leakage Other Symptoms: \_\_\_\_\_

4. Detail of Request  Repair (Overhaul)  Regular Checks

5. Others: \_\_\_\_\_

Company Name: \_\_\_\_\_ Personnel in charge: \_\_\_\_\_

Address: \_\_\_\_\_

Tel: \_\_\_\_\_ Fax: \_\_\_\_\_ E-mail: \_\_\_\_\_

Agent Name: \_\_\_\_\_ Personnel in charge: \_\_\_\_\_

Address: \_\_\_\_\_

Tel: \_\_\_\_\_ Fax: \_\_\_\_\_

\* In case you do not have any direct transaction with us, please be sure to fill in the agent name.

6. Confirmation

The gas and substance used in this pump or unit is harmless to human bodies, or it is not contaminated by any substance harmful to human bodies.

Signed \_\_\_\_\_ (seal) Date: \_\_\_\_\_

\* In order to avoid a trouble during transportation, please evacuate oil from any oil pump before shipping.

\* You are requested to ship the package to our Service Division (CS Center). (See the attached list of addresses.)

# Sales, service agency, and the where to make contact

## <Head Office/ Miyazaki Plan>

291-7 Chausubaru, Saito-shi, Miyazaki Prf. Japan  
TEL (81)983-42-1411 FAX (81) 983-42-1422

## <Sales Office>

Sales Division & Overseas Division & Yokohama Branch  
1-10-4, Kitashinyokohama, Kohoku-ku, Yokohama-shi, Kanagawa Prf. Japan

Yokohama branch: TEL (81) 45-533-0203 FAX (81) 45-533-0204

Overseas Division: TEL (81) 45-533-0206 FAX (81) 45-533-0204

## Osaka Branch

3-3-31 Miyahara, Yodogawa-ku, Osaka-shi, Osaka-fu Japan  
TEL(81)6-6350-2166 FAX(81)6-6350-2169

## Nagoya Branch

3-11-31 Sakae, Naka-ku, Nagoya-shi, Aichi Prf. Japan  
TEL(81)52-249-5121 FAX(81)52-249-5122

## <Service Office>

### Yokohama Branch

1-10-4, Kitashinyokohama, Kohoku-ku, Yokohama-shi, Kanagawa Prf. Japan  
TEL(81)45-533-0509 FAX(81)45-533-0512

### Miyazaki Branch

291-7 Chausubaru, Saito-shi, Miyazaki Prf. Japan  
TEL(81)983-42-4135 FAX(81)983-43-2159

## ULVAC GmbH

Parkring 11, 85748, Garching ,Germany  
TEL(49)89-96-0909-0 FAX(49)89-96-0909-96